



# SZABO SCANDIC

Part of Europa Biosite

## Produktinformation



Forschungsprodukte & Biochemikalien



Zellkultur & Verbrauchsmaterial



Diagnostik & molekulare Diagnostik



Laborgeräte & Service

Weitere Information auf den folgenden Seiten!  
See the following pages for more information!



### Lieferung & Zahlungsart

siehe unsere [Liefer- und Versandbedingungen](#)

### Zuschläge

- Mindermengenzuschlag
- Trockeneiszuschlag
- Gefahrgutzuschlag
- Expressversand

### SZABO-SCANDIC HandelsgmbH

Quellenstraße 110, A-1100 Wien

T. +43(0)1 489 3961-0

F. +43(0)1 489 3961-7

[mail@szabo-scandic.com](mailto:mail@szabo-scandic.com)

[www.szabo-scandic.com](http://www.szabo-scandic.com)

[linkedin.com/company/szaboscandic](https://www.linkedin.com/company/szaboscandic) 

## **Data Sheet**

### ***TDO Inhibitor Screening Assay Kit (384)***

**Catalog # 72036**

**BACKGROUND:** L-tryptophan (L-Trp) is an essential amino acid necessary for protein synthesis in mammalian cells and the L-Trp to kynurenine (Kyn) pathway is firmly established as a key regulator of innate and adaptive immunity. Catabolism of L-Trp to Kyn maintains an immunosuppressive microenvironment by starving immune cells of L-Trp and releasing degradation products of L-Trp that have immunosuppressive functions. Tryptophan 2,3 dioxygenase (TDO) is upregulated in many tumors, providing cancer cells with an avenue for immune evasion.

**DESCRIPTION:** The *TDO Inhibitor Screening Assay Kit* is designed to measure TDO enzyme inhibition. The kit comes in a convenient format, with enough reaction solution and enzyme to perform a total of 400 reactions. The *TDO Inhibitor Screening Assay Kit* is simple to use. Inhibitor and enzyme are added to a sample containing L-Trp substrate. After a room temperature incubation, activity is measured by measuring the absorption of reaction product at  $\lambda = 320\text{-}325\text{ nm}$ .

**COMPONENTS:**

Catalog #	Component	Amount	Storage	
71195	TDO, His-Tag	100 $\mu\text{g}$	-80°C	<b>(Avoid freeze/thaw cycles!)</b>
73005	TDO Reaction Solution	4 x 10 ml	-80°C	
73006	TDO Assay Buffer	2 x 1 ml	-80°C	
	UV transparent 384-well plate	1	Room Temp.	

**MATERIALS OR INSTRUMENTS REQUIRED BUT NOT SUPPLIED:**

Spectrophotometer capable of measuring absorbance at  $\lambda = 320\text{-}325\text{ nm}$  in a 384 well plate.

**APPLICATIONS:** Useful for the study of TDO enzymology, screening inhibitors and selectivity profiling.

**CONTRAINDICATIONS:** DMSO >1%, strong acids or bases, ionic detergents, high salt

**STABILITY:** Stable at least 6 months from date of receipt, when stored as directed. Kit components require different storage conditions. Be sure to store each component at the proper temperature upon arrival.

**REFERENCE:** Li, J.S., *et al. Arch. Insect Biochem. Physiol.* 2007 Feb; **64(2)**: 74–87.

OUR PRODUCTS ARE FOR RESEARCH USE ONLY. NOT FOR DIAGNOSTIC OR THERAPEUTIC USE.

To place your order, please contact us by Phone **1.858.202.1401**, Fax **1.858.481.8694**  
Or you can Email us at: [info@bpsbioscience.com](mailto:info@bpsbioscience.com)  
Please visit our website at: [www.bpsbioscience.com](http://www.bpsbioscience.com)

**ASSAY PROTOCOL:**

All samples and controls should be tested in duplicate. Use slow shaking for all incubations.

**Step 1:**

- 1) Thaw **TDO Reaction Solution** and aliquot 90  $\mu$ l into each well. *Note: **TDO Reaction Solution** may contain a precipitate after thawing. Please ensure the mixture is fully solubilized by gentle mixing before aliquoting. Do not vortex.*
- 2) Add 5  $\mu$ l of inhibitor solution (no more than 20% DMSO) to each well designated "Test Inhibitor". For the "Positive Control" and "Blank", add 5  $\mu$ l of the same solution without inhibitor (inhibitor buffer). *Note: Keep the final DMSO concentration below 1%.*
- 3) Thaw **TDO enzyme** on ice. Upon first thaw, briefly spin tube containing enzyme to recover full contents of the tube. Aliquot **TDO enzyme** into single use aliquots. Store remaining undiluted enzyme in aliquots at -80°C. *Note: **TDO enzyme** is very sensitive to freeze/thaw cycles. Do not re-use thawed aliquots or diluted enzyme.*
- 4) Dilute **TDO** in **TDO Assay Buffer** at 50 ng/ $\mu$ l. Keep diluted protein on ice until use. Discard any unused diluted protein after use.



	Blank	Positive Control	Test Inhibitor
Reaction Solution	90 $\mu$ l	90 $\mu$ l	90 $\mu$ l
Test Inhibitor	–	–	5 $\mu$ l
Inhibitor buffer (no inhibitor)	5 $\mu$ l	5 $\mu$ l	–
TDO Assay Buffer	5 $\mu$ l	–	–
TDO (50 ng/ $\mu$ l)	–	5 $\mu$ l	5 $\mu$ l
<b>Total</b>	<b>100 <math>\mu</math>l</b>	<b>100 <math>\mu</math>l</b>	<b>100 <math>\mu</math>l</b>

- 5) Add 5  $\mu$ l of **TDO Assay Buffer** to the well designated "Blank".
- 6) Initiate reaction by adding 5  $\mu$ l of diluted **TDO** prepared as described above to the wells labeled "Positive Control", and "Test Inhibitor". Incubate at room temperature for 90 minutes.
- 7) Measure absorption at  $\lambda$ =320-325 nm. Subtract "Blank" value from all other values.

OUR PRODUCTS ARE FOR RESEARCH USE ONLY. NOT FOR DIAGNOSTIC OR THERAPEUTIC USE.

To place your order, please contact us by Phone **1.858.202.1401**, Fax **1.858.481.8694**  
Or you can Email us at: [info@bpsbioscience.com](mailto:info@bpsbioscience.com)  
Please visit our website at: [www.bpsbioscience.com](http://www.bpsbioscience.com)

**EXAMPLE OF ASSAY RESULTS:**

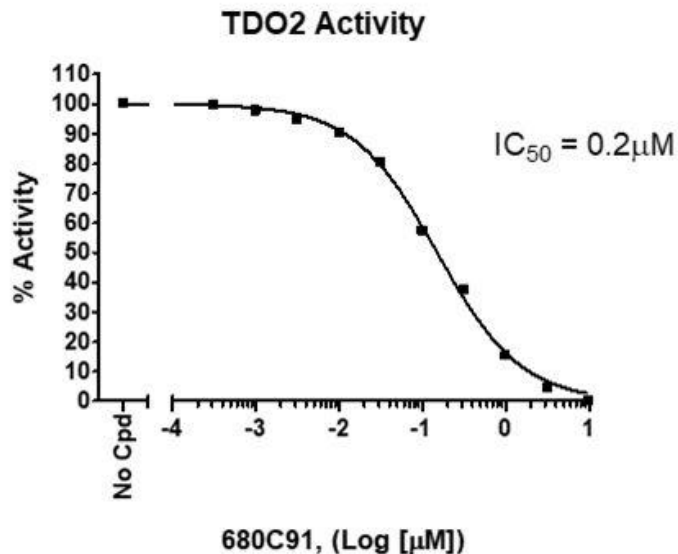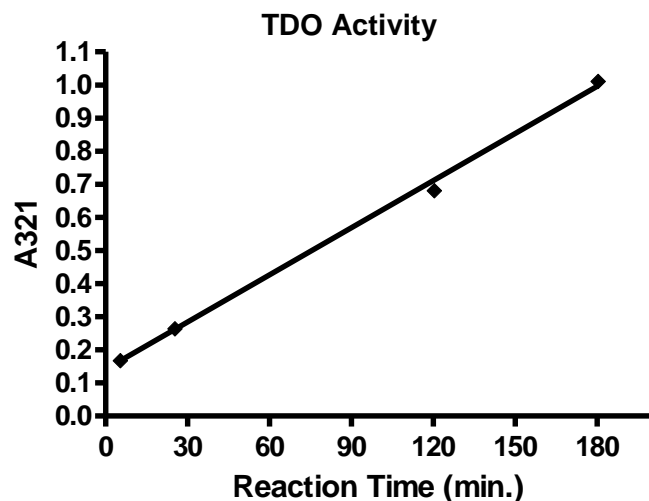

TDO activity (left) and inhibition (right), measured using the TDO Inhibitor Screening Assay Kit, BPS Bioscience, Catalog #73026. Data shown is lot-specific. For lot-specific information, please contact BPS Bioscience, Inc. at [info@bpsbioscience.com](mailto:info@bpsbioscience.com).

OUR PRODUCTS ARE FOR RESEARCH USE ONLY. NOT FOR DIAGNOSTIC OR THERAPEUTIC USE.

To place your order, please contact us by Phone **1.858.202.1401**, Fax **1.858.481.8694**  
 Or you can Email us at: [info@bpsbioscience.com](mailto:info@bpsbioscience.com)  
 Please visit our website at: [www.bpsbioscience.com](http://www.bpsbioscience.com)



6042 Cornerstone Ct. West, Ste. B  
San Diego, CA 92121  
**Tel:** 1.858.202.1401  
**Fax:** 1.858.481.8694  
**Email:** [info@bpsbioscience.com](mailto:info@bpsbioscience.com)

**RELATED PRODUCTS:**

<b><u>Product Name</u></b>	<b><u>Catalog#</u></b>	<b><u>Size</u></b>
TDO, His-tag	71195	50 µg
IDO1, His-tag	71182	50 µg
IDO2, His-tag	71194	50 µg
IDO1 Inhibitor Screening Assay Kit	72021	96 rxns
IDO1 Inhibitor Screening Assay Kit	72034	384 rxns
IDO2 Inhibitor Screening Assay Kit	72022	96 rxns
TDO Inhibitor Screening Assay Kit	72023	96 rxns
IDO1 Fluorogenic Inhibitor Screening Assay Kit	72037	96 rxns
TDO Fluorogenic Inhibitor Screening Assay Kit	72039	96 rxns
IDO1 Cell-Based Assay Kit	72031	100 rxns
TDO Cell-Based Assay Kit	72033	100 rxns
NLG919	27337-1	10 mg
INCB024360	27338-1	10 mg
IDO1 Reaction Solution	73001	10 ml
IDO1 Assay Buffer	73002	1 ml
IDO2 Reaction Solution	73003	10 ml
IDO2 Assay Buffer	73004	1 ml
TDO Reaction Solution	73005	10 ml
TDO Assay Buffer	73006	1 ml

OUR PRODUCTS ARE FOR RESEARCH USE ONLY. NOT FOR DIAGNOSTIC OR THERAPEUTIC USE.

To place your order, please contact us by Phone **1.858.202.1401**, Fax **1.858.481.8694**  
Or you can Email us at: [info@bpsbioscience.com](mailto:info@bpsbioscience.com)  
Please visit our website at: [www.bpsbioscience.com](http://www.bpsbioscience.com)