



SZABO SCANDIC

Part of Europa Biosite

Produktinformation



Forschungsprodukte & Biochemikalien



Zellkultur & Verbrauchsmaterial



Diagnostik & molekulare Diagnostik



Laborgeräte & Service

Weitere Information auf den folgenden Seiten!
See the following pages for more information!



Lieferung & Zahlungsart

siehe unsere [Liefer- und Versandbedingungen](#)

Zuschläge

- Mindermengenzuschlag
- Trockeneiszuschlag
- Gefahrgutzuschlag
- Expressversand

SZABO-SCANDIC HandelsgmbH

Quellenstraße 110, A-1100 Wien

T. +43(0)1 489 3961-0

F. +43(0)1 489 3961-7

mail@szabo-scandic.com

www.szabo-scandic.com

[linkedin.com/company/szaboscandic](https://www.linkedin.com/company/szaboscandic) 

Description

The STAT3 Luciferase Reporter THP-1 cell line is designed for monitoring the STAT3 signal transduction pathway. It contains a firefly luciferase gene driven by STAT3 response elements located upstream of the minimal TATA promoter. After activation by cytokines or growth factors, endogenous STAT3 binds to the DNA response elements, inducing transcription of the luciferase reporter gene.

Background

The JAK-STAT signaling pathway is involved in both cancer and autoimmune disease and has been an attractive target for drug discovery. In the context of immune regulation, STAT3 is expressed in diverse cell types. While it was initially described as an acute phase response factor in the context of IL-6 signaling, STAT3 has since been identified as a downstream regulator of many different cytokines. STAT3-mediated IL-6 signaling in macrophages has also been shown to play a role in their differentiation and proliferation. Studies of the JAK-STAT signaling pathways can prove useful in furthering our understanding of immunological diseases and cancer.

Application

- Monitor the STAT3 signaling pathway in a monocyte cell line
- Screen for compound activity on the STAT3 signaling pathway
- Screen for inhibitors of STAT3-mediated IL-6 signaling

Materials Provided

| Components | Format |
|-------------------------|-------------------------------------------------------------------------------------------------------|
| 2 vials of frozen cells | Each vial contains $\sim 2 \times 10^6$ cells in 1 ml of cell freezing medium (BPS Bioscience #79796) |

Parental Cell Line

THP-1, monocyte, suspension

Mycoplasma Testing

The cell line has been screened to confirm the absence of Mycoplasma species.

Materials Required but Not Supplied

These materials are not supplied with the cell line but are necessary for cell culture and cellular assays. BPS Bioscience's reagents are validated and optimized for use with this cell line and are highly recommended for best results. Media components are provided in the Media Formulations section below.

Media Required for Cell Culture

| Name | Ordering Information |
|------------------|---------------------------------------|
| Thaw Medium 8 | BPS Bioscience #79652 |
| Growth Medium 8A | BPS Bioscience #79653 |

Materials Required for Cellular Assay

| Name | Ordering Information |
|-----------------------------------|---------------------------------------|
| Thaw Medium 8 | BPS Bioscience #79652 |
| Recombinant Human IL-6 Protein | R&D Systems #206-IL |
| Pan-JAK inhibitor CP 690,550 | Cayman #11598 |
| ONE-Step™ Luciferase Assay System | BPS Bioscience #60690 |
| Luminometer | |

Storage Conditions



Cells are shipped in dry ice and should immediately be thawed or stored in liquid nitrogen upon receipt. Do not use a -80°C freezer for long term storage. Contact technical support at support@bpsbioscience.com if the cells are not frozen in dry ice upon arrival.

Media Formulations

For best results, it is *highly recommended* to use these validated and optimized media from BPS Bioscience. Other preparations or formulations of media may result in suboptimal performance.



Note: Thaw Media do *not* contain selective antibiotics. However, Growth Media *do* contain selective antibiotics, which are used for maintaining the presence of the transfected gene(s) over passages. Cells should be grown at 37°C with 5% CO₂. BPS Bioscience's cell lines are stable for at least 15 passages when grown under proper conditions.

Media Required for Cell Culture

Thaw Medium 8 (BPS Bioscience #79652):

RPMI 1640 medium supplemented with 10% heat-inactivated FBS, 1% Penicillin/Streptomycin

Growth Medium 8A (BPS Bioscience #79653):

RPMI 1640 medium supplemented with 10% heat-inactivated FBS, 1% Penicillin/Streptomycin plus 1 µg/ml of Puromycin.

Media Required for Functional Cellular Assay

Thaw Medium 8 (BPS Bioscience #79652):

RPMI 1640 medium supplemented with 10% heat-inactivated FBS, 1% Penicillin/Streptomycin

Cell Culture Protocol

Cell Thawing

1. Swirl the vial of frozen cells for approximately 60 seconds in a 37°C water bath. As soon as the cells are thawed (it may be slightly faster or slower than 60 seconds), quickly transfer the entire contents of the vial to a tube containing 10 ml of pre-warmed Thaw Medium 8 (**no Puromycin**).

Leaving the cells in the water bath at 37°C for too long will result in rapid loss of viability.

2. Immediately spin down the cells at 300 x g for 5 minutes, remove the medium and resuspend the cells in 5 ml of pre-warmed Thaw Medium 8 (**no Puromycin**).
3. Transfer the resuspended cells to a T25 flask and incubate at 37°C in a 5% CO₂ incubator.
4. After 24 hours of culture, check for cell viability. For a T25 flask, add 3-4 ml of Thaw Medium 8 (**no Puromycin**), and continue growing in a 5% CO₂ incubator at 37°C until the cells are ready to passage.
5. Cells should be passaged before they reach a density of 2 x 10⁶ cells/ml. At first passage and subsequent passages, use Growth Medium 8A (**contains Puromycin**).

Cell Passage

Dilute the cell suspension into new culture vessels before they reach a density of 2 x 10⁶ cells/ml, at no less than 0.2 x 10⁶ cells/ml of Growth Medium 8A (**contains Puromycin**). The sub-cultivation ratio should maintain the cells between 0.5 x 10⁶ cells/ml and 2 x 10⁶ cells/ml.

Cell Freezing

1. Spin down the cells at 300 x g for 5 minutes, remove the medium and resuspend the cell pellet in 4°C Freezing Medium (BPS Bioscience #79796, or 10% DMSO + 90% FBS) at a density of $\sim 2 \times 10^6$ cells/ml.
2. Dispense 1 ml of cell aliquots into cryogenic vials. Place the vials in an insulated container for slow cooling and store at -80°C overnight.
3. Transfer the vials to liquid nitrogen the next day for storage.



Note: It is recommended to expand the cells and freeze at least 10 vials at an early passage for future use.

Assay Performance

Functional characterization of STAT3 Luciferase Reporter THP-1 Cell Line.

The following assays are designed for 96-well format. To perform the assay in different tissue culture formats, the cell number and reagent volumes should be scaled appropriately.

Assay Medium:

Thaw Medium 8 (BPS Bioscience #79652):

RPMI 1640 medium supplemented with 10% heat-inactivated FBS, 1% Penicillin/Streptomycin

A. Human IL-6 dose response

1. Seed STAT3 Luciferase Reporter THP-1 cells at a density of 60,000 cells per well into a white, clear-bottom 96-well cell culture plate in 90 μ l of Thaw medium 8. Incubate the cells at 37°C with 5% CO₂ overnight.
2. Set up each treatment at least in triplicates:

Prepare a 3-fold increment serial dilution of human IL-6 in Thaw medium 8 at 10X the desired final concentrations (prepare enough for 10 μ l/well).

- a. Add 10 μ l of each dilution to wells labeled "hIL-6 stimulated wells".
 - b. Add 10 μ l of Thaw medium 8 to unstimulated control wells (for measuring uninduced level of STAT3 reporter activity)
 - c. Add 100 μ l of Thaw medium 8 to cell-free control wells (for determining background luminescence).
3. Incubate at 37°C with 5% CO₂ for \sim 5-6 hours
 4. Perform luciferase assay using ONE-Step™ Luciferase Assay System according to the protocol provided:
 - a. Add 100 μ l of ONE-Step™ Luciferase reagent per well and rock at room temperature for \sim 15 minutes.

- b. Measure luminescence using a luminometer. If using luciferase reagents from other vendors, follow the manufacturer's assay protocol.
5. Data Analysis: Subtract the average background luminescence (cell-free control wells) from the luminescence reading of all conditions. The fold induction of STAT3 luciferase reporter expression is the background-subtracted luminescence of hIL-6-stimulated cells divided by the background-subtracted luminescence of unstimulated control cells.

$$\text{Fold induction} = \frac{(\text{luminescence of stimulated cells} - \text{background})}{(\text{luminescence of unstimulated cells} - \text{background})}$$

Validation Data

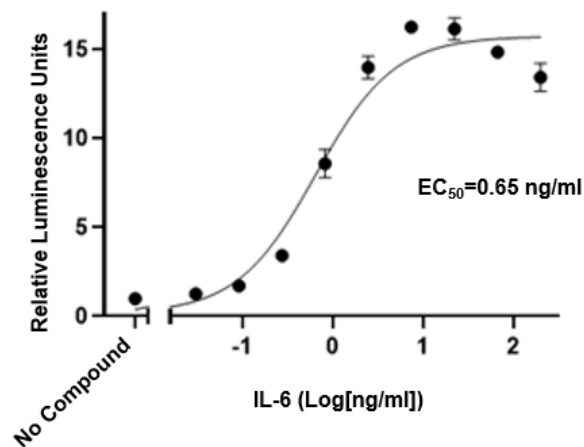


Figure 1: Dose response of human IL-6 in STAT3 Luciferase Reporter THP-1 cells.

Increasing concentrations of human IL-6 were added to the cells the day after seeding. Controls consisted of untreated cells and no-cell wells (background). Approximately 5 hours later, the ONE-Step™ luciferase assay reagent was added to the cells. The results are shown as fold induction of luciferase reporter expression. Fold induction was determined by comparing values of IL-6-stimulated cells against the value for control cells without hIL-6 treatment. The EC₅₀ of hIL-6 in this cell line was 0.65 ng/ml.

B. Inhibition of IL-6-induced STAT3 Reporter Activity by JAK inhibitor CP 690,550 in STAT3 Luciferase Reporter THP-1 cells

1. Seed the STAT3 Luciferase Reporter THP-1 cells at a density of 60,000 cells per well into white, clear-bottom 96-well cell culture plate in 90 µl of Thaw medium 8.
2. Set up each treatment in at least triplicates:

Prepare a 3-fold increment serial dilution of CP 690,550 in Thaw medium 8 at 10X the desired concentrations (prepare enough for 10 µl/well).

- a. Add 10 µl of each dilution to wells labeled "CP 690,550 treated wells".

- b. Add 10 μ l of Thaw medium 8 to untreated control wells (for measuring uninduced level of STAT3 reporter activity)
3. Incubate the cells at 37°C with 5% CO₂ overnight.
4. Stimulate the cells with hIL-6:
 - a. Add 10 μ l of hIL-6 diluted in Thaw medium 8 to the cells except for control unstimulated cells. The recommended final hIL-6 concentration is 10 ng/ml.
 - b. Add 10 μ l of Thaw medium 8 to the unstimulated control cells (for determining STAT3 basal activity).
 - c. Add 110 μ l of Thaw medium 8 to cell-free control wells (for determining background luminescence).
5. Incubate at 37°C with 5% CO₂ for ~5-6 hours
6. Perform luciferase assay using ONE-Step™ Luciferase Assay System according to the protocol provided.
7. Add 110 μ l of ONE-Step™ Luciferase reagent per well and rock at room temperature for ~15 minutes.
8. Measure luminescence using a luminometer. If using other luciferase reagents from other vendors, follow the manufacturer's assay protocol.
9. Data Analysis: Subtract the average background luminescence (cell-free control wells) from the average luminescence reading of all conditions. The Relative Luminescence value is the background-subtracted luminescence of CP 690,550-treated cells divided by the background-subtracted luminescence of untreated cells, reported to 100.

$$\text{Relative luminescence} = \left(\frac{\text{luminescence of drug treated cells} - \text{background}}{\text{luminescence of untreated cells} - \text{background}} \right) \times 100$$

Validation Data

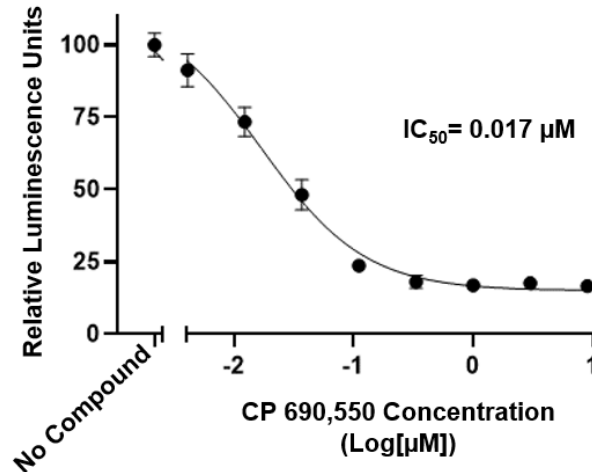


Figure 2: Inhibition of hIL-6-induced STAT3 Reporter Activity by JAK inhibitor CP 690,550 in STAT3 Luciferase Reporter THP-1 cells.

The cells were incubated with increasing concentrations of pan-JAK inhibitor CP 690,550 overnight, then stimulated with human IL-6 (10 ng/ml) for approximately 5 hours. Controls consisted of untreated cells (no compound) and no-cell wells (background). Luciferase activity was measured using the ONE-Step™ Luciferase assay system. The results are shown as percentage of luminescence of STAT3 reporter activity (in which IL-6-stimulated cells in the absence of JAK inhibitor is set at 100%). The IC₅₀ of CP 690,550 in this cell line was 0.017 μM.

References

- Akira S., *et al* (1994). *Cell* **77**(1):63-71
 Hammaren H. M., *et al.* (2019). *Cytokine* **118**:48-63
 Hillmer E., *et al.* (2016). *Cytokine & Growth Factor Review* **31**: 1-15
 Nakajima K., *et al.*, (1996). *EMBO J.* **14**: 3651-8
 Niemand C., *et al.* (2003). *J. Immunology.* **170**(6): 3263-3272

License Disclosure

Visit bpsbioscience.com/license for the label license and other key information about this product.

Troubleshooting Guide

Visit bpsbioscience.com/cell-line-faq for detailed troubleshooting instructions. For all further questions, please email support@bpsbioscience.com.

Related Products

| <i>Products</i> | <i>Catalog #</i> | <i>Size</i> |
|-----------------------------------------------------|------------------|---------------|
| STAT3 Reporter Jurkat Cell Line | 78497 | 500 µl x 2 |
| STAT3 eGFP Reporter Lentivirus | 78197 | 500 µl x 2 |
| STAT3 Luciferase Reporter Lentivirus | 79744 | 500 µl x 2 |
| STAT3 Reporter (Luc) - HEK293 Cell line (Puromycin) | 79800-P | 2 vials |
| STAT3 Reporter Kit (STAT3 Signaling Pathway) | 79730 | 500 reactions |
| STAT3, GST-tag Recombinant | 75003 | 20 µg |
| Human Interferon-alpha 2a Recombinant | 90158-A | 20 µg |