



SZABO SCANDIC

Part of Europa Biosite

Produktinformation



Forschungsprodukte & Biochemikalien



Zellkultur & Verbrauchsmaterial



Diagnostik & molekulare Diagnostik



Laborgeräte & Service

Weitere Information auf den folgenden Seiten!
See the following pages for more information!



Lieferung & Zahlungsart

siehe unsere [Liefer- und Versandbedingungen](#)

Zuschläge

- Mindermengenzuschlag
- Trockeneiszuschlag
- Gefahrgutzuschlag
- Expressversand

SZABO-SCANDIC HandelsgmbH

Quellenstraße 110, A-1100 Wien

T. +43(0)1 489 3961-0

F. +43(0)1 489 3961-7

mail@szabo-scandic.com

www.szabo-scandic.com

[linkedin.com/company/szaboscandic](https://www.linkedin.com/company/szaboscandic) 



6042 Cornerstone Court W, Ste B
San Diego, CA 92121
Tel: 1.858.829.3082
Fax: 1.858.481.8694
Email: info@bpsbioscience.com

Data Sheet
ABL (T315I) Assay Kit
Catalog #79576
96 reactions

DESCRIPTION: The Abelson (ABL) family of nonreceptor tyrosine kinases is critical to the progression and treatment of chronic myeloid leukemia (CML). The single-point mutation T315I has been linked to drug resistance and poor prognosis, making it an important drug target. The *ABL (T315I) Assay Kit* is designed to measure ABL activity for screening and profiling applications using Kinase-Glo[®] MAX as a detection reagent. The *ABL (T315I) Assay Kit* comes in a convenient 96-well format, with enough purified recombinant ABL enzyme, ABLtide peptide, ATP, and kinase assay buffer for 100 enzyme reactions.

COMPONENTS:

Catalog #	Reagent	Amount	Storage	
40415	ABL (T315I)	1 µg	-80°C	Avoid multiple freeze/thaw cycles!
	5x Kinase assay buffer	1.5 ml	-20°C	
	ATP (500 µM)	100 µl	-20°C	
79577	ABLtide (1 mg/ml)	500 µl	-20°C	
	96-well plate, white	1	Room Temp.	

MATERIALS OR INSTRUMENTS REQUIRED BUT NOT SUPPLIED:

Kinase-Glo MAX (Promega #V6071)
Dithiothreitol (DTT, 1 M; optional)
Microplate reader capable of reading luminescence
Adjustable micropipettor and sterile tips
30°C incubator

APPLICATIONS: Useful for studying enzyme kinetics and screening small molecular inhibitors for drug discovery and HTS applications.

STABILITY: Up to 6 months when stored as recommended.

REFERENCE:

1. Greuber, E. K., Smith-Pearson, P., Wang, J., & Pendergast, A. M. (2013). Role of ABL Family Kinases in Cancer: from Leukemia to Solid Tumors. *Nature Reviews. Cancer*, **13(8)**, 559–571.
2. Ganguly SS, Plattner R. (2012). Activation of Abl Family Kinases in Solid Tumors. *Genes & Cancer* **3(5-6)**:414-425.

OUR PRODUCTS ARE FOR RESEARCH USE ONLY. NOT FOR DIAGNOSTIC OR THERAPEUTIC USE.

To place your order, please contact us by Phone **1.858.829.3082** Fax **1.858.481.8694**

Or you can Email us at: info@bpsbioscience.com

Please visit our website at: www.bpsbioscience.com



6042 Cornerstone Court W, Ste B
San Diego, CA 92121
Tel: 1.858.829.3082
Fax: 1.858.481.8694
Email: info@bpsbioscience.com

ASSAY PROTOCOL:

All samples and controls should be tested in duplicate.

- 1) Thaw **5x Kinase assay buffer**, **ATP (500 μ M)**, and **ABLtide (1 mg/ml)**.
(Optional: If desired, add DTT to **5x Kinase assay buffer** to make a 10 mM concentration; e.g. add 10 μ l of 1 M DTT to 1 ml **5x Kinase assay buffer**)
- 2) Prepare the master mixture (25 μ l per well): N wells x (5 μ l **5x Kinase assay buffer** + 1 μ l **ATP (500 μ M)** + 5 μ l **ABLtide (1 mg/ml)** + 14 μ l water). Add 25 μ l to every well.

	Positive Control	Test Inhibitor	Blank
5x Kinase assay buffer	5 μ l	5 μ l	5 μ l
ATP (500 μ M)	1 μ l	1 μ l	1 μ l
ABLtide(1 mg/ml)	5 μ l	5 μ l	5 μ l
Water	14 μ l	14 μ l	14 μ l
Test Inhibitor	–	5 μ l	–
Inhibitor Buffer (no inhibitor)	5 μ l	–	5 μ l
1x Kinase buffer	–	–	20 μ l
ABL (T315I) (0.5 ng/ μ l)	20 μ l	20 μ l	–
Total	50 μ l	50 μ l	50 μ l

- 3) Add 5 μ l of Inhibitor solution of each well labeled as "Test Inhibitor." For the "Positive Control" and "Blank," add 5 μ l of the same solution without inhibitor (Inhibitor buffer). *Note: Keep DMSO concentration of the Test Inhibitor at $\leq 10\%$, as final DMSO concentration in the reaction should be $\leq 1\%$.*
- 4) Prepare 3 ml of **1x Kinase assay buffer** by mixing 600 μ l of **5x Kinase assay buffer** with 2400 μ l water. 3 ml of **1x Kinase assay buffer** is sufficient for 100 reactions.
- 5) To the wells designated as "Blank," add 20 μ l of **1x Kinase assay buffer**.
- 6) Thaw **ABL (T315I)** on ice. Upon first thaw, briefly spin tube containing enzyme to recover full content of the tube. Calculate the amount of **ABL (T315I)** required for the assay and dilute enzyme to 0.5 ng/ μ l with **1x Kinase assay buffer**. Store remaining undiluted enzyme in aliquots at -80°C . *Note: ABL (T315I) enzyme is sensitive to freeze/thaw cycles. Avoid multiple freeze/thaw cycles. Do not re-use thawed aliquots or diluted enzyme.*

OUR PRODUCTS ARE FOR RESEARCH USE ONLY. NOT FOR DIAGNOSTIC OR THERAPEUTIC USE.

To place your order, please contact us by Phone **1.858.829.3082** Fax **1.858.481.8694**

Or you can Email us at: info@bpsbioscience.com

Please visit our website at: www.bpsbioscience.com



6042 Cornerstone Court W, Ste B
San Diego, CA 92121
Tel: 1.858.829.3082
Fax: 1.858.481.8694
Email: info@bpsbioscience.com

- 7) Initiate reaction by adding 20 μ l of diluted **ABL (T315I)** to the wells designated "Positive Control" and "Test Inhibitor Control." Incubate at 30°C for 40 minutes.
- 8) Thaw Kinase-Glo Max reagent.
- 9) After the 40 minute reaction, add 50 μ l of Kinase-Glo Max reagent to each well. Cover plate with aluminum foil and incubate the plate at room temperature for 15 minutes.
- 10) Immediately read sample in a luminometer or microtiter-plate capable of reading chemiluminescence. "Blank" value is subtracted from all readings.

Reading Chemiluminescence:

Chemiluminescence is the emission of light (luminescence) which results from a chemical reaction. The detection of chemiluminescence requires no wavelength selection because the method used is emission photometry and is not emission spectrophotometry.

To properly read chemiluminescence, make sure the plate reader is set for LUMINESCENCE mode. Typical integration time is 1 second, delay after plate movement is 100 msec. Do not use a filter when measuring light emission. Typical settings for the Synergy 2 BioTek plate reader are: use the "hole" position on the filter wheel; Optics position: Top; Read type: endpoint. Sensitivity may be adjusted based on the luminescence of a control assay without enzyme (typically we set this value as 100).

OUR PRODUCTS ARE FOR RESEARCH USE ONLY. NOT FOR DIAGNOSTIC OR THERAPEUTIC USE.

To place your order, please contact us by Phone **1.858.829.3082** Fax **1.858.481.8694**


Or you can Email us at: info@bpsbioscience.com

Please visit our website at: www.bpsbioscience.com



6042 Cornerstone Court W, Ste B
San Diego, CA 92121
Tel: 1.858.829.3082
Fax: 1.858.481.8694
Email: info@bpsbioscience.com

Example of Assay Results:

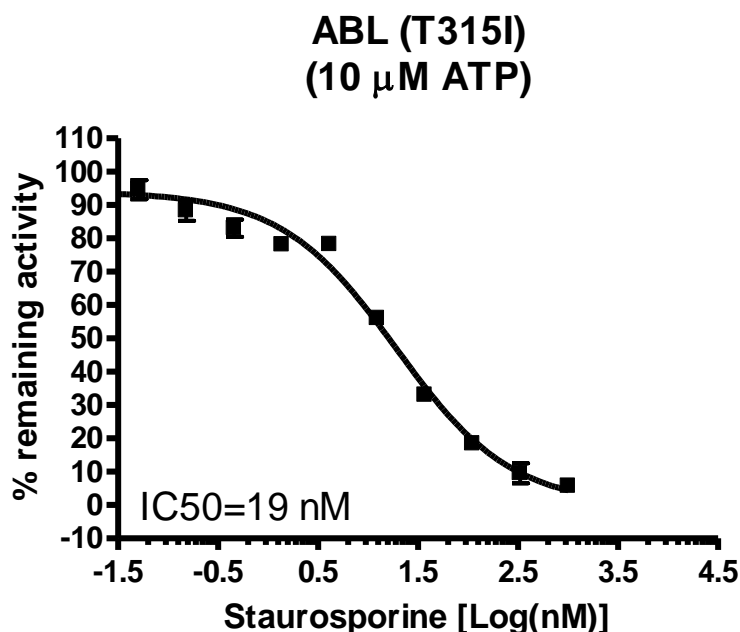

Inhibition of ABL enzyme by Staurosporine (BPS Bioscience, #27002), measured using the ABL (T315I) assay kit (Cat. #79576). *Data shown is lot-specific. For lot-specific information, please contact BPS Bioscience, Inc. at info@bpsbioscience.com*

RELATED PRODUCTS:

<u>Product Name</u>	<u>Catalog #</u>	<u>Size</u>
Abl Kinase, His-tag	40411	10 μ g
Abl Kinase Mutant (T315I), His-tag	40415	10 μ g
Abl2, His-tag	40417	10 μ g
Staurosporine	27002	10 mg
5x Kinase assay buffer	79334	10 ml

OUR PRODUCTS ARE FOR RESEARCH USE ONLY. NOT FOR DIAGNOSTIC OR THERAPEUTIC USE.

To place your order, please contact us by Phone **1.858.829.3082** Fax **1.858.481.8694**

Or you can Email us at: info@bpsbioscience.com

Please visit our website at: www.bpsbioscience.com