



SZABO SCANDIC

Part of Europa Biosite

Produktinformation



Forschungsprodukte & Biochemikalien



Zellkultur & Verbrauchsmaterial



Diagnostik & molekulare Diagnostik



Laborgeräte & Service

Weitere Information auf den folgenden Seiten!
See the following pages for more information!



Lieferung & Zahlungsart

siehe unsere [Liefer- und Versandbedingungen](#)

Zuschläge

- Mindermengenzuschlag
- Trockeneiszuschlag
- Gefahrgutzuschlag
- Expressversand

SZABO-SCANDIC HandelsgmbH

Quellenstraße 110, A-1100 Wien

T. +43(0)1 489 3961-0

F. +43(0)1 489 3961-7

mail@szabo-scandic.com

www.szabo-scandic.com

[linkedin.com/company/szaboscandic](https://www.linkedin.com/company/szaboscandic)



DCAF1 Intrachain TR-FRET Assay Kit

Description

The DCAF1 Intrachain TR-FRET Assay Kit is a homogeneous, sensitive TR-FRET (Time-Resolved Fluorescence Resonance Energy Transfer) assay kit designed to measure DCAF1 (DDB1 and CUL4-associated Factor 1) complex auto-ubiquitination activity. It utilizes a mix of donor and acceptor labeled ubiquitin to complete the TR-FRET pairing. This assay measures poly-ubiquitination since both the TR-FRET donor and acceptor are incorporated into poly-ubiquitin chains formed on DCAF1. The kit contains enough purified DCAF1/Rbx1 (M5I)/DDB1/CUL4A Complex, purified UBE1 and UbcH5b, detection reagents and assay buffer for 400 reactions.

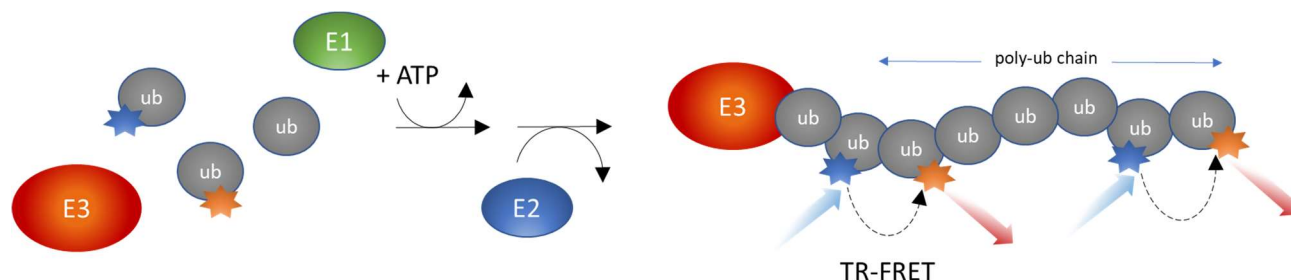


Figure 1: DCAF1 Intrachain TR-FRET Assay Kit schematic.

The TR-FRET signal is proportional to DCAF1 auto-polyubiquitination activity.

Background

Covalent conjugation to ubiquitin (Ub) is one of the major post-translational modifications that regulates protein stability, function, and localization. Ubiquitination is the concerted action of three enzymes: a Ub-activating enzyme (E1), a Ub-conjugating enzyme (E2), and a Ub ligase (E3). The specificity and efficiency of ubiquitination are largely determined by the E3 enzyme, which directs the last step of the Ub-conjugating cascade by binding to both an E2~Ub conjugate and a substrate protein. This step ensures the transfer of Ub from E2~Ub to the substrate, leading to its mono- or poly-ubiquitination.

DCAF1 (DDB1 and CUL4 Associated Factor 1), also known as VprBP (Vpr binding protein), is a WD40 repeat domain-containing E3 ligase of the CRL4 (cullin RING ligase 4) subfamily of proteins. It functions as the substrate recognition unit in E3 ligase complexes, such as the EED (embryonic ectoderm development)-DDB1 (DNA damage-binding protein)-VprBP (or EDVP) and the CRL4^{DCAF1} complex. The CRL4^{DCAF1} complex is composed of CUL4A (cullin-4A), DDB1 and DCAF1. DCAF1 is involved in various normal physiological functions in the cell, including cell cycle regulation/progression, cell division, lipid metabolism, and miRNA biogenesis. Additionally, DCAF1 plays an E3 ligase-independent role, by acting as a kinase and phosphorylating Histone H2A. Antagonists of DCAF1 could be used toward the development of therapeutics for cancers and viral treatments. DCAF1 is found at high levels in cancer cells, and it is implicated in disease progression. The use of a DCAF1-BTK (Bruton's tyrosine kinase) PROTAC[®] resulted in degradation of BTK in cells that had acquired resistance to CRBN (cereblon)-BTK PROTACs, making DCAF1 an interesting alternative E3 ligase in cases of resistance mechanism to other commonly used PROTACs. A deeper understanding of the role of this protein and how to manipulate its multiple activities may open new therapeutic avenues.

Applications

- Screen molecules that inhibit DCAF1 E3 ligase activity in drug discovery High-Throughput Screening (HTS) applications.
- Determine the IC₅₀ of inhibitors of DCAF1 E3 ligase activity.
- Determine DCAF1 E3 ligase activity real-time kinetics.

Supplied Materials

Catalog #	Name	Amount	Storage
80301	UBE1, FLAG-Tag*	50 µg	-80°C
80314	UbcH5b, His-Tag (Human)*	60 µg	-80°C
102190	DCAF1/Rbx1 (M5I)/DDB1/CUL4A Complex*	32 µg	-80°C
82185	200x Ubi-Mix™	40 µl	-80°C
	4 mM ATP	2 x 1 ml	-80°C
78856	U2 Assay Buffer	2 x 10 ml	-80°C
79969	White, nonbinding, low volume 384-well microtiter plate	1	Room Temp

* The initial concentration of enzyme is lot-specific and will be indicated on the tube containing the protein.

Materials Required but Not Supplied

- Fluorescent microplate reader capable of measuring Time Resolved Fluorescence Resonance Energy Transfer (TR-FRET)
- Adjustable micropipettor and sterile tips
- Orbital Shaker

Storage Conditions

This assay kit will perform optimally for up to **6 months** from date of receipt when the materials are stored as directed.

Safety

This product is for research purposes only and not for human or therapeutic use. This product should be considered hazardous and is harmful by inhalation, in contact with skin, eyes, clothing, and if swallowed. If contact occurs, wash thoroughly.

Contraindications

- This kit is compatible with up to 1% final DMSO concentration.
- Compounds that are fluorescent may interfere with the results, depending on their spectral excitation and emission properties.
- It is recommended that the compound alone is tested to determine any potential interference of the compound on the assay results.

Assay Protocol

- All samples and controls should be performed in triplicate.
- The assay should include “Blank”, “Positive Control”, “Negative Control” and “Test Inhibitor” conditions.
- We recommend maintaining the diluted protein on ice during use.
- For detailed information on protein handling please refer to Protein FAQs (bpsbioscience.com).

- We recommend using Methylated Ubiquitin Recombinant (#102075) as internal control. If not running a dose response curve for the control inhibitor, we recommend running the control inhibitor at 0.1X, 1X and 10X the IC₅₀ value shown in the validation data below.
- For instructions on how to prepare reagent dilutions please refer to [Serial Dilution Protocol \(bpsbioscience.com\)](https://bpsbioscience.com).

1. Thaw **UBE1**, **Ubch5b**, **DCAF1 complex**, **U2 Assay Buffer**, **200 x Ubi-Mix™**, and **ATP** on ice. Briefly spin the tubes to recover their full content.
2. Dilute 200x Ubi-Mix™ 40-fold with U2 Assay Buffer. This makes 5x Ubi-Mix™.

Note: Do not reuse the diluted solution. 200 x Ubi-Mix™ is sensitive to freeze/thaw cycles. Prepare single use aliquots (minimum volume of 5 µl/aliquot) and freeze immediately at -80°C.

3. Dilute **UBE1** to 96 ng/µl with U2 Assay Buffer (1 µl/well).
4. Dilute **Ubch5b** to 144 ng/µl with U2 Assay Buffer (1 µl/well).
5. Dilute **DCAF1 complex** to 16 ng/µl with U2 Assay Buffer (5 µl/well).
6. Prepare the Test Inhibitor (4 µl/well): for a titration, prepare serial dilutions at concentrations 5-fold higher than the desired final concentrations. The final volume of the reaction is 20 µl.

6.1 If the Test Inhibitor is water-soluble, prepare serial dilutions in U2 Assay Buffer, 5-fold more concentrated than the desired final concentrations.

For the positive and negative controls, use U2 Assay Buffer (Diluent Solution).

OR

6.2 If the Test inhibitor is soluble in DMSO, prepare the test inhibitor at a concentration 100-fold higher than the highest desired concentration in 100% DMSO, then dilute the inhibitor 20-fold in U2 Assay Buffer to prepare the highest concentration of the 5-fold intermediate dilutions. The concentration of DMSO is now 5%.

Prepare serial dilutions of the Test Inhibitor at 5-fold the desired final concentrations using 5% DMSO in U2 Assay Buffer to keep the concentration of DMSO constant.

For controls prepare 5% DMSO in U2 Assay Buffer (Diluent Solution) so that all wells contain the same amount of DMSO.

Note: The final concentration of DMSO should not exceed 1%.

7. Prepare a Deficient Master Mix (15 µl/ "Blank" well): N wells x (4 µl of 5x Ubi-Mix™ + 1 µl of diluted UBE1 + 1 µl of diluted Ubch5b + 4 µl of Diluent Solution + 5 µl of Assay Buffer).

8. Add 15 µl of Deficient Master Mix to the “Blank” wells.
9. Prepare a Master Mix (11 µl/well, except “Blank” wells): N wells x (4 µl of 5x Ubi-Mix™ + 1 µl of diluted UBE1 + 1 µl of diluted Ubch5b + 5 µl of diluted DCAF1 complex).
10. Add 11 µl of Master Mix to all wells, except the “Blank” wells.
11. Add 4 µl of diluted inhibitor to the “Test Inhibitor” wells.
12. Add 4 µl of Diluent Solution to the “Positive Control” and “Negative Control” wells.
13. Initiate the reaction by adding 5 µl of **4 mM ATP** to the “Blank”, “Test Inhibitor”, and “Positive Control” wells.
14. Add 5 µl of U2 Assay Buffer to the “Negative Control” wells.
15. Read the fluorescent intensity in a microtiter-plate reader capable of measuring TR-FRET in kinetic mode for up to 1 hour. An end point readout can be done in 30-40 minutes.

	Blank	Test Inhibitor	Positive Control	Negative Control
Deficient Master Mix	15 µl	-	-	-
Master Mix	-	11 µl	11 µl	11 µl
U2 Assay Buffer	-	-	-	5 µl
Test Inhibitor	-	4 µl	-	-
Diluent Solution	-	-	4 µl	4 µl
4 mM ATP	5 µl	5 µl	5 µl	-
Total	20 µl	20 µl	20 µl	20 µl

Instrument Settings

Two sequential measurements should be conducted. Eu- donor emission should be measured at 620 nm followed by dye-acceptor emission at 665 nm.

Reading Mode	Time Resolved
Excitation Wavelength	317 (20 nm bandwidth)
Emission Wavelength	620 (10 nm bandwidth)
Lag Time	60 µs
Integration Time	500 µs
Excitation Wavelength	317 (20 nm bandwidth)
Emission Wavelength	665 (10 nm bandwidth)
Lag Time	60 µs
Integration Time	500 µs

CALCULATING RESULTS

Data analysis is performed using the TR-FRET ratio (665 nm emission/620 nm emission).

$$FRET = \frac{S_{665}}{S_{620}}$$

When percentage activity is calculated, the FRET value from the Blank (it is expected that Blank and Negative Control have a similar values) can be set as zero percent activity and the FRET value from the positive control can be set as one hundred percent activity.

$$\% \text{ Activity} = \frac{FRET_S - FRET_{blank}}{FRET_P - FRET_{blank}} \times 100\%$$

FRET_S = FRET value for samples of Test Inhibitor, FRET_{blank} = FRET value for the Blank, and FRET_P = FRET value for the Positive Control (no inhibitor).

Example Results

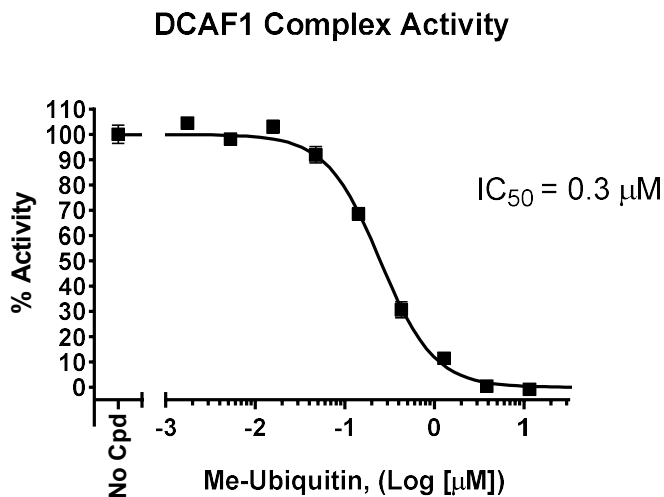


Figure 2: Inhibition of DCAF1 auto-ubiquitination activity by Methylated Ubiquitin Recombinant. DCAF1 auto-ubiquitination was measured in presence of increasing concentrations of Methylated Ubiquitin Recombinant (#102075). Results are expressed as percent activity, in which absence of inhibitor is set to 100%.

Data shown is representative. For lot-specific information, please contact BPS Bioscience, Inc. at support@bpsbioscience.com.

Troubleshooting Guide

Visit bpsbioscience.com/assay-kits-faq for detailed troubleshooting instructions. For all further questions, please email support@bpsbioscience.com

References

Li A., *et al.*, 2023 *J Med Chem* 66 (7): 5041-5060.

Schröder M., *et al.*, 2024 *Nat Commun* 15 (1): 275.

Related Products

<i>Products</i>	<i>Catalog #</i>	<i>Size</i>
Cereblon Intrachain TR-FRET Assay Kit	78301	384 reactions
MDM2 Intrachain TR-FRET Assay Kit	78302	384 reactions
SMURF1 Intrachain TR-FRET Assay Kit	78303	384 reactions
VHL Intrachain TR-FRET Assay Kit	78305	384 reactions
DCAF15 Intrachain TR-FRET Assay Kit	78543	384 reactions
DCAF11 Intrachain TR-FRET Assay Kit	78542	384 reactions

Version 052124