



SZABO SCANDIC

Part of Europa Biosite

Produktinformation



Forschungsprodukte & Biochemikalien



Zellkultur & Verbrauchsmaterial



Diagnostik & molekulare Diagnostik



Laborgeräte & Service

Weitere Information auf den folgenden Seiten!
See the following pages for more information!



Lieferung & Zahlungsart

siehe unsere [Liefer- und Versandbedingungen](#)

Zuschläge

- Mindermengenzuschlag
- Trockeneiszuschlag
- Gefahrgutzuschlag
- Expressversand

SZABO-SCANDIC HandelsgmbH

Quellenstraße 110, A-1100 Wien

T. +43(0)1 489 3961-0

F. +43(0)1 489 3961-7

mail@szabo-scandic.com

www.szabo-scandic.com

[linkedin.com/company/szaboscandic](https://www.linkedin.com/company/szaboscandic) 

Description

The Chemi-Verse™ HER2 Kinase Assay Kit is designed to measure HER2 (human epidermal growth factor receptor 2) tyrosine kinase activity for screening and profiling applications using ADP-Glo™ as a detection reagent. The assay kit comes in a convenient 96-well format, with enough purified HER2 (amino acids 679-1255), kinase substrate, ATP, and kinase assay buffer for 100 enzyme reactions.

Background

HER2 (human epidermal growth factor receptor 2), also known as erbB-2 or CD340, is a tyrosine kinase of the EGFR family of proteins. There is no known ligand, but it can form homodimers or heterodimers with other HER proteins. Once active, it activates the MAPK (mitogen-activated protein kinase) and PI3K (phosphatidylinositol-3 kinase) signaling pathways resulting in cell cycle progression and cell proliferation. HER2 over-expression is also known to occur in breast, ovarian, stomach, lung adenocarcinoma, aggressive forms of uterine cancer and gastric cancer. In 1990 the FDA approved the use of the monoclonal antibody trastuzumab in breast and stomach cancer. Other strategies to target HER2 that have been approved include ADCs (antibody-drug conjugate) and margetuximab (an HER2 antibody that can alter the Fc-receptor affinity to CD16 and induce cytotoxicity). The use of small molecule tyrosine kinase inhibitors, alone or in combinatory therapy, has shown great promise in the treatment of HER2⁺ breast cancer (BC). However, resistance to treatment, for instance by mutations on HER2 or upregulation of other HER receptors, has been described. Neratinib, a pan-HER2 inhibitor, was approved in 2017 for early-stage BC, as adjuvant anti-HER2 therapy after trastuzumab treatment. However, side effects limit its use. The development of treatments able to target early-stage cancer, with minimal side effects and resistance development, will bring major benefits to HER2⁺ oncology patients.

Applications

Study enzyme kinetics and screen small molecule inhibitors for drug discovery and high throughput screening (HTS) applications.

Supplied Materials

Catalog #	Name	Amount	Storage
40230	HER2, GST-Tag*	15 µg	-80°C
79334	5x Kinase Buffer 1	1.5 ml	-20°C
79686	500 µM ATP	50 µl	-20°C
40217	PTK substrate Poly(Glu:Tyr 4:1) (10 mg/ml)	50 µl	-20°C
82545	White 96-well plate	1	Room Temperature

*The concentration of the protein is lot-specific and will be indicated on the tube.

Materials Required but Not Supplied

Name	Ordering Information
ADP-Glo™ Kinase Assay	Promega #V6930
DTT (Dithiothreitol), 1M, optional	
Microplate reader capable of reading luminescence	
Adjustable micropipettor and sterile tips	
30°C incubator	

Storage Conditions

This assay kit will perform optimally for up to **6 months** from date of receipt when the materials are stored as directed.

Safety

This product is for research purposes only and not for human or therapeutic use. This product should be considered hazardous and is harmful by inhalation, in contact with skin, eyes, clothing, and if swallowed. If contact occurs, wash thoroughly.

Assay Principle

The **ADP-Glo™ Kinase Assay (Promega #V6930)** quantifies the amount of ADP produced by a kinase upon phosphorylation of a substrate. First, addition of the ADP-Glo™ reagent terminates the reaction and quenches the remaining ATP. Second, the addition of the Kinase Detection reagent converts the produced ADP to ATP. The newly generated ATP is quantified by a luciferase reaction. The luminescent signal correlates with the amount of ADP generated by the kinase and is linear to 1 mM ATP.

Contraindications

The final concentration of DMSO in the assay should not exceed 1%.

Assay Protocol

- All samples and controls should be tested in duplicate.
- The assay should include “Blank”, “Positive Control” and “Test Inhibitor” conditions.
- We recommend maintaining the diluted protein on ice during use.
- For detailed information on protein handling please refer to [Protein FAQs \(bpsbioscience.com\)](https://www.bpsbioscience.com/protein-faqs).
- We recommend using Afatinib (#27009) or Epertinib hydrochloride (#82559) as internal control. If not running a dose response curve for the control inhibitor, we recommend running the control inhibitor at 0.1X, 1X and 10X the IC₅₀ value shown in the validation data below.
- For instructions on how to prepare reagent dilutions please refer to [Serial Dilution Protocol \(bpsbioscience.com\)](https://www.bpsbioscience.com/serial-dilution-protocol).

1. Thaw **5x Kinase Assay Buffer 1**, **500 μM ATP**, and **PTK substrate Poly(Glu:Tyr 4:1)(10 mg/ml)**.

Optional: If desired, make 5x Kinase Assay Buffer 1 with 10 mM DTT.

2. Prepare 3 ml of **1x Kinase Assay Buffer 1** by mixing 600 μl of **5x Kinase Assay Buffer 1** with 2,400 μl of distilled water.

Note: Three (3 ml) of 1x Kinase Assay Buffer 1 is sufficient for 100 reactions.

3. Prepare a **Master Mix** (12.5 μl/well): N wells x (6 μl of 5x Kinase Assay Buffer 1 + 0.5 μl of 500 μM ATP + 0.5 μl of PTK substrate Poly(Glu:Tyr 4:1)(10 mg/ml) + 5.5 μl of distilled water).
4. Add 12.5 μl of Master Mix to every well.

5. Prepare the **Test Inhibitor** (2.5 µl/well): for a titration prepare serial dilutions at concentrations 10-fold higher than the desired final concentrations. The final volume of the reaction is 25 µl.

5.1 If the Test Inhibitor is water-soluble: Prepare serial dilutions in 1x Kinase Assay Buffer 1, 10-fold more concentrated than the desired final concentrations.

For the positive and negative controls, use 1x Kinase Assay Buffer 1 (Diluent Solution).

OR

5.2 If the Test inhibitor is soluble in DMSO: Prepare the test inhibitor at 100-fold the highest desired concentration in 100% DMSO, then dilute the inhibitor 10-fold in 1x Kinase Assay Buffer 1 to prepare the highest concentration of the 10-fold intermediate dilutions. The concentration of DMSO is now 10%.

Prepare serial dilutions of the Test Inhibitor at 10-fold the desired final concentrations using 10% DMSO in 1x Kinase Assay Buffer 1 to keep the concentration of DMSO constant.

For positive and negative controls, prepare 10% DMSO in 1x Kinase Assay Buffer 1 (vol/vol) so that all wells contain the same amount of DMSO (Diluent Solution).

Note: The final concentration of DMSO should not exceed 1%.

6. Add 2.5 µl of Test Inhibitor to each well labeled "Test Inhibitor".
7. Add 2.5 µl of Diluent Solution to the "Positive Control" and "Blank" wells.
8. Add 10 µl of 1x Kinase Assay Buffer 1 to the "Blank" wells.
9. Thaw **HER2 Kinase** on ice. Briefly spin the tube to recover its full content.
10. Dilute the protein kinase (10 µl/well) to 14 ng/µl with **1x Kinase Assay Buffer 1**.
11. Initiate the reaction by adding 10 µl of diluted kinase to the wells designated "Positive Control" and "Test Inhibitor".

Component	Blank	Positive Control	Test Inhibitor
Master Mix	12.5 µl	12.5 µl	12.5 µl
Test Inhibitor	-	-	2.5 µl
Diluent Solution	2.5 µl	2.5 µl	-
1x Kinase Assay Buffer 1	10 µl	-	-
Diluted HER2 (14 ng/µl)	-	10 µl	10 µl
Total	25 µl	25 µl	25 µl

12. Incubate at 30°C for 45 minutes.

13. Thaw the ADP-Glo™ reagent.
14. At the end of the 45 minute reaction, add 25 µl of ADP-Glo™ reagent to each well.
15. Cover the plate with aluminum foil and incubate at Room Temperature (RT) for 45 minutes.
16. Thaw the Kinase Detection Reagent.
17. Add 50 µl of Kinase Detection reagent to each well.
18. Cover the plate with aluminum foil and incubate at RT for another 45 minutes.
19. Immediately read in a luminometer or a microplate reader capable of reading luminescence.
20. The “Blank” value is subtracted from all other readings.

Reading Luminescence

Luminescence is the emission of light resulting from a chemical reaction. The detection of luminescence requires no wavelength selection because the method used is emission photometry and not emission spectrophotometry.

To properly read luminescence, make sure the plate reader is set for LUMINESCENCE mode. Typical integration time is 1 second, delay after plate movement is 100 msec. Do not use a filter when measuring light emission. Typical settings for the Synergy 2 BioTek plate reader: use the “hole” position on the filter wheel; Optics position: Top; Read type: endpoint. Sensitivity may be adjusted based on the luminescence of a control assay without enzyme (typically we set this value as 100).

Example Results

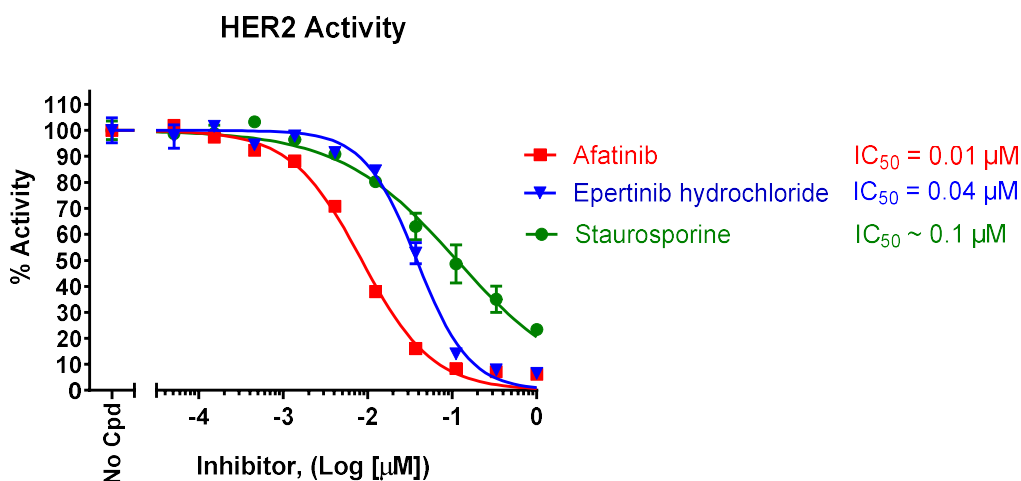


Figure 1: Inhibition of HER2 kinase activity by the inhibitors Afatinib, Epertinib hydrochloride and Staurosporine.

HER2 kinase activity was measured in the presence of increasing concentrations of Afatinib (#27009), Epertinib hydrochloride (#82559) or Staurosporine (#27002). The “Blank” value was subtracted from all other values. Results are expressed as the percent of control (kinase activity in the absence of inhibitor, set at 100%).

Data shown is representative. For lot-specific information, please contact BPS Bioscience, Inc. at support@bpsbioscience.com.

Troubleshooting Guide

Visit bpsbioscience.com/assay-kits-faq for detailed troubleshooting instructions. For all further questions, please email support@bpsbioscience.com

References

Schlam I. and Swain S., 2021 *npj Breast Cancer* 7: 56.

Related Products

Products	Catalog #	Size
Chemi-Verse™ HER4 Kinase Assay Kit	82553	96 reactions
HER2 (ERBB2) CHO Recombinant Cell Line (High, Medium, or Low Expression)	79612	2 Vials
Afatinib	27009	25 mg
Staurosporine	27002	1 mg
HER3 CHO Cell Line	82299	2 vials
HER3, FC Fusion, Avi-Tag Recombinant	102228	20 μg/50 μg

Version 053124