



SZABO SCANDIC

Part of Europa Biosite

Produktinformation



Forschungsprodukte & Biochemikalien



Zellkultur & Verbrauchsmaterial



Diagnostik & molekulare Diagnostik



Laborgeräte & Service

Weitere Information auf den folgenden Seiten!
See the following pages for more information!



Lieferung & Zahlungsart

siehe unsere [Liefer- und Versandbedingungen](#)

Zuschläge

- Mindermengenzuschlag
- Trockeneiszuschlag
- Gefahrgutzuschlag
- Expressversand

SZABO-SCANDIC HandelsgmbH

Quellenstraße 110, A-1100 Wien

T. +43(0)1 489 3961-0

F. +43(0)1 489 3961-7

mail@szabo-scandic.com

www.szabo-scandic.com

[linkedin.com/company/szaboscandic](https://www.linkedin.com/company/szaboscandic) 



Progesterone ELISA Kit

Item No. 582601

www.caymanchem.com

Customer Service 800.364.9897

Technical Support 888.526.5351

1180 E. Ellsworth Rd • Ann Arbor, MI • USA

TABLE OF CONTENTS

GENERAL INFORMATION	3	Materials Supplied
	4	Safety Data
	4	Precautions
	5	If You Have Problems
	5	Storage and Stability
	5	Materials Needed but Not Supplied
INTRODUCTION	6	Background
	6	About This Assay
	7	Description of AChE Competitive ELISAs
	8	Biochemistry of Acetylcholinesterase
	10	Definition of Key Terms
PRE-ASSAY PREPARATION	11	Buffer Preparation
	12	Sample Preparation
	13	Sample Purification
	15	Sample Matrix Properties
ASSAY PROTOCOL	17	Preparation of Assay-Specific Reagents
	19	Plate Set Up
	20	Performing the Assay
ANALYSIS	23	Calculations
	26	Performance Characteristics
RESOURCES	30	Troubleshooting
	31	Additional Reading
	31	References
	33	Plate Template
	34	Notes
	35	Warranty and Limitation of Remedy

GENERAL INFORMATION

Materials Supplied

Item Number	Item	96 wells Quantity/Size	480 wells Quantity/Size
482602	Progesterone ELISA Antiserum	1 vial/100 dtn	1 vial/500 dtn
482600	Progesterone AChE Tracer	1 vial/100 dtn	1 vial/500 dtn
482604	Progesterone ELISA Standard	1 vial	1 vial
400060	ELISA Buffer Concentrate (10X)	2 vials/10 ml	4 vials/10 ml
400062	Wash Buffer Concentrate (400X)	1 vial/5 ml	1 vial/12.5 ml
400035	Polysorbate 20	1 vial/3 ml	1 vial/3 ml
400004/400006	Mouse Anti-Rabbit IgG Coated Plate	1 plate	5 plates
400012	96-Well Cover Sheet	1 cover	5 covers
400050	Ellman's Reagent	3 vials/100 dtn	6 vials/250 dtn

If any of the items listed above are damaged or missing, please contact our Customer Service department at (800) 364-9897 or (734) 971-3335. We cannot accept any returns without prior authorization.



WARNING: THIS PRODUCT IS FOR RESEARCH ONLY - NOT FOR HUMAN OR VETERINARY DIAGNOSTIC OR THERAPEUTIC USE.

Safety Data

This material should be considered hazardous until further information becomes available. Do not ingest, inhale, get in eyes, on skin, or on clothing. Wash thoroughly after handling. Before use, the user must review the complete Safety Data Sheet, which has been sent *via* email to your institution.

Precautions

Please read these instructions carefully before beginning this assay.

The reagents in this kit have been tested and formulated to work exclusively with Cayman's AChE ELISA Kits. This kit may not perform as described if any reagent or procedure is replaced or modified.

When compared to quantification by LC/MS or GC/MS, it is not uncommon for immunoassays to report higher analyte concentrations. While LC/MS or GC/MS analyses typically measure only a single compound, antibodies used in immunoassays sometimes recognize not only the target molecule, but also structurally related molecules, including biologically relevant metabolites. In many cases, measurement of both the parent molecule and metabolites is more representative of the overall biological response than is the measurement of a short-lived parent molecule. It is the responsibility of the researcher to understand the limits of both assay systems and to interpret their data accordingly.

If You Have Problems

Technical Service Contact Information

Phone: 888-526-5351 (USA and Canada only) or 734-975-3888
Fax: 734-971-3640
Email: techserv@caymanchem.com

In order for our staff to assist you quickly and efficiently, please be ready to supply the lot number of the kit (found on the outside of the box).

Storage and Stability

This kit will perform as specified if stored as directed at -20°C and used before the expiration date indicated on the outside of the box.

Materials Needed But Not Supplied

1. A plate reader capable of measuring absorbance between 405-420 nm.
2. An adjustable pipettor.
3. A source of 'UltraPure' water. Water used to prepare all ELISA reagents and buffers must be deionized and free of trace organic contaminants ('UltraPure'). Use activated carbon filter cartridges or other organic scavengers. Glass distilled water (even if double distilled), HPLC-grade water, and sterile water (for injections) are not adequate for ELISA. *NOTE: UltraPure water is available for purchase from Cayman (Item No. 400000).*
4. Materials used for Sample Preparation (see page 12).

Background

Progesterone, along with pregnenolone, is the biosynthetic precursor of all other steroid hormones. Progesterone is synthesized from cholesterol by the sequential action of desmolase in the mitochondria, which produces pregnenolone, followed by $\Delta^{4,5}$ -isomerase in the outer mitochondrial membrane and smooth endoplasmic reticulum of steroid-secreting cells.¹ The main function of progesterone is to prepare the uterine lining for implantation of a fertilized ovum and to maintain pregnancy. Measurement of serum or plasma progesterone levels is used as an index to monitor ovulation and investigate luteal function. Plasma concentrations of progesterone are approximately 0.2-0.8 ng/ml and 4-20 ng/ml during the follicular phase and luteal phase of the menstrual cycle, respectively.¹ Salivary concentrations of progesterone are approximately 100-fold lower than those found in plasma.²

About This Assay

Cayman's Progesterone ELISA Kit is a competitive assay that can be used for quantification of progesterone in plasma and other sample matrices. The assay has a range from 7.8-1,000 pg/ml and a sensitivity (80% B/B₀) of approximately 10 pg/ml.

Description of AChE Competitive ELISAs^{3,4}

This assay is based on the competition between progesterone and a progesterone-acetylcholinesterase (AChE) conjugate (progesterone tracer) for a limited number of progesterone-specific rabbit antiserum binding sites. Because the concentration of the progesterone tracer is held constant while the concentration of progesterone varies, the amount of progesterone tracer that is able to bind to the rabbit antiserum will be inversely proportional to the concentration of progesterone in the well. This rabbit antiserum-progesterone (either free or tracer) complex binds to the mouse monoclonal anti-rabbit IgG that has been previously attached to the well. The plate is washed to remove any unbound reagents and then Ellman's Reagent (which contains the substrate to AChE) is added to the well. The product of this enzymatic reaction has a distinct yellow color and absorbs strongly at 412 nm. The intensity of this color, determined spectrophotometrically, is proportional to the amount of progesterone tracer bound to the well, which is inversely proportional to the amount of free progesterone present in the well during the incubation; or

$$\text{Absorbance} \propto [\text{Bound Progesterone Tracer}] \propto 1/[\text{Progesterone}]$$

A schematic of this process is shown in Figure 1, below.

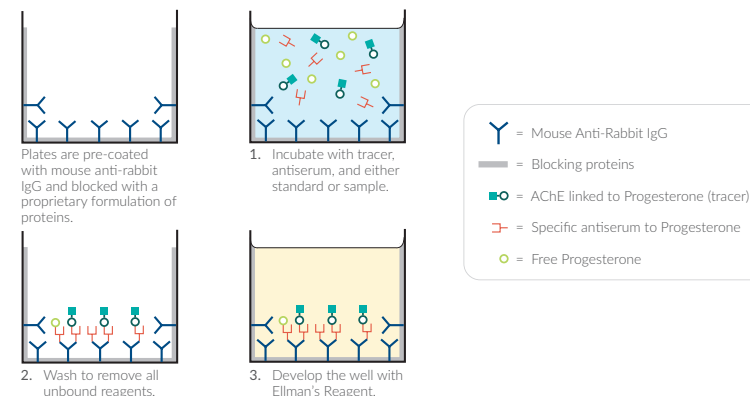


Figure 1. Schematic of the AChE ELISA

Biochemistry of Acetylcholinesterase

The electric organ of the electric eel, *E. electricus*, contains an avid AChE capable of massive catalytic turnover during the generation of its electrochemical discharges. The electric eel AChE has a clover leaf-shaped tertiary structure consisting of a triad of tetramers attached to a collagen-like structural fibril. This stable enzyme is capable of high turnover ($64,000 \text{ s}^{-1}$) for the hydrolysis of acetylthiocholine.

A molecule of the analyte covalently attached to a molecule of AChE serves as the tracer in AChE enzyme immunoassays. Quantification of the tracer is achieved by measuring its AChE activity with Ellman's Reagent. This reagent consists of acetylthiocholine and 5,5'-dithio-bis-(2-nitrobenzoic acid). Hydrolysis of acetylthiocholine by AChE produces thiocholine (see Figure 2, page 9). The non-enzymatic reaction of thiocholine with 5,5'-dithio-bis-(2-nitrobenzoic acid) produces 5-thio-2-nitrobenzoic acid, which has a strong absorbance at 412 nm ($\epsilon = 13,600$).

AChE has several advantages over other enzymes commonly used for enzyme immunoassays. Unlike horseradish peroxidase, AChE does not self-inactivate during turnover. This property of AChE also allows redevelopment of the assay if it is accidentally splashed or spilled. In addition, the enzyme is highly stable under the assay conditions, has a wide pH range (pH 5-10), and is not inhibited by common buffer salts or preservatives. Since AChE is stable during the development step, it is unnecessary to use a "stop" reagent, and the plate may be read whenever it is convenient.

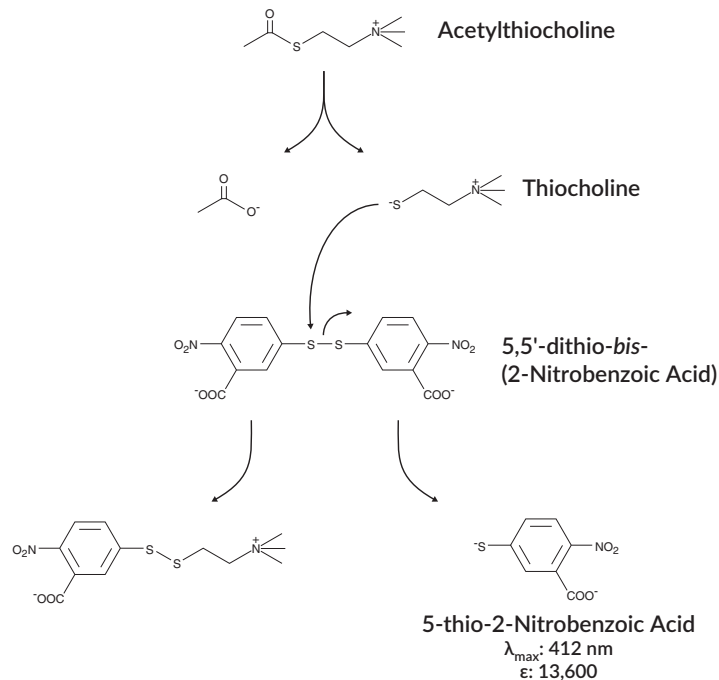


Figure 2. Reaction catalyzed by acetylcholinesterase

Definition of Key Terms

Blank: background absorbance caused by Ellman's Reagent. The blank absorbance should be subtracted from the absorbance readings of all the other wells, including NSB wells.

Total Activity: total enzymatic activity of the AChE-linked tracer. This is analogous to the specific activity of a radioactive tracer.

NSB (Non-Specific Binding): non-immunological binding of the tracer to the well. Even in the absence of specific antibody a very small amount of tracer still binds to the well; the NSB is a measure of this low binding. Do not forget to subtract the Blank absorbance values.

B₀ (Maximum Binding): maximum amount of the tracer that the antibody can bind in the absence of free analyte.

%B/B₀ (%Bound/Maximum Bound): ratio of the absorbance of a particular sample or standard well to that of the maximum binding (B₀) well.

Standard Curve: a plot of the %B/B₀ values *versus* concentration of a series of wells containing various known amounts of analyte.

Dtn: determination, where one dtn is the amount of reagent used per well.

Cross Reactivity: numerical representation of the relative reactivity of this assay towards structurally related molecules as compared to the primary analyte of interest. Biomolecules that possess similar epitopes to the analyte can compete with the assay tracer for binding to the primary antibody. Substances that are superior to the analyte in displacing the tracer result in a cross reactivity that is greater than 100%. Substances that are inferior to the primary analyte in displacing the tracer result in a cross reactivity that is less than 100%. Cross reactivity is calculated by comparing the mid-point (50% B/B₀) value of the tested molecule to the mid-point (50% B/B₀) value of the primary analyte when each is measured in assay buffer using the following formula:

$$\% \text{ Cross Reactivity} = \left[\frac{50\% \text{ B/B}_0 \text{ value for the primary analyte}}{50\% \text{ B/B}_0 \text{ value for the potential cross reactant}} \right] \times 100\%$$

PRE-ASSAY PREPARATION

NOTE: Water used to prepare all ELISA reagents and buffers must be deionized and free of trace organic contaminants ('UltraPure'). Use activated carbon filter cartridges or other organic scavengers. Glass distilled water (even if double distilled), HPLC-grade water, and sterile water (for injections) are not adequate for ELISA. UltraPure water may be purchased from Cayman (Item No. 400000).

Buffer Preparation

Store all diluted buffers at 4°C; they will be stable for about two months.

1. ELISA Buffer Preparation

Dilute the contents of one vial of ELISA Buffer Concentrate (10X) (Item No. 400060) with 90 ml of UltraPure water. Be certain to rinse the vial to remove any salts that may have precipitated. *NOTE: It is normal for the concentrated buffer to contain crystalline salts after thawing. These will completely dissolve upon dilution with water.*

2. Wash Buffer Preparation

5 ml vial Wash Buffer Concentrate (400X) (96-well kit; Item No. 400062): Dilute to a total volume of 2 liters with UltraPure water and add 1 ml of Polysorbate 20 (Item No. 400035).

OR

12.5 ml vial Wash Buffer Concentrate (400X) (480-well kit; Item No. 400062): Dilute to a total volume of 5 liters with UltraPure water and add 2.5 ml of Polysorbate 20 (Item No. 400035).

Smaller volumes of Wash Buffer can be prepared by diluting the Wash Buffer Concentrate 1:400 and adding Polysorbate 20 (0.5 ml/liter of Wash Buffer).

NOTE: Polysorbate 20 is a viscous liquid and cannot be measured by a regular pipette. A positive displacement pipette or a syringe should be used to deliver small quantities accurately.

Sample Preparation

General Precautions

- All samples must be free of organic solvents prior to assay.
- Samples should be assayed immediately after collection; samples that cannot be assayed immediately should be stored at -80°C.
- Samples of rabbit origin may contain antibodies which interfere with the assay by binding to the mouse anti-rabbit plate. We recommend that all rabbit samples be purified prior to use in this assay.
- AEBSF (Pefabloc SC®) and PMSF inhibit AChE. Samples containing these protease inhibitors should not be used in this assay.
- Fetal bovine serum (FBS) may contain significant amounts of steroids that could interfere with this assay. We recommend using FBS that has been charcoal-stripped or dialyzed to remove free hormone.

Sample Purification

In general, tissue culture supernatant samples may be diluted with ELISA Buffer and added directly to the assay well.

Testing for Interference

Plasma and serum, as well as other heterogeneous mixtures such as lavage fluids and aspirates often contain contaminants which can interfere in the assay. It is best to check for interference before embarking on a large number of sample measurements. To test for interference, dilute one or two test samples to obtain at least two different dilutions of each sample between 10 and 500 pg/ml (i.e. between ~20-80% B/B₀). If the two dilutions of the sample show good correlation (differ by 20% or less) in the final calculated progesterone concentration, purification is not required. If you do not see good correlation of the different dilutions, purification is advised. The **Purification Protocol** on page 14 is one such method.

NOTE: We do not recommend using heparin in plasma collection for this assay.

Plasma and Serum

The following protocol is a suggestion only. You may choose a different protocol based on your own requirements, sample type and expertise.

1. Aliquot a known amount of each sample into a clean test tube (500 µl is recommended). If your samples need to be concentrated, a larger volume should be used (e.g., a 5 ml sample will be concentrated by a factor of 10, a 10 ml sample will be concentrated by a factor of 20, etc.).
2. Add 4X the sample volume of 3% isopropanol in hexane, or Methyl tert-butyl ether (MTBE), or diethyl ether (DEE), and mix thoroughly with a vortexer. Allow layers to separate. Transfer the organic (upper) layer into a clean tube using a transfer pipette. Repeat this extraction procedure two to five times.*
3. Evaporate the organic solvent under a gentle stream of nitrogen.
4. To resuspend the sample, add 500 µl of ELISA Buffer. Vortex. It is common for insoluble precipitate to remain in the sample after the addition of ELISA Buffer; this will not affect the assay. This sample is now ready for use in the ELISA.

*If it is necessary to stop during this purification, samples may be stored in the organic solvent at -20°C or -80°C.

Sample Matrix Properties

Spike and Recovery

Plasma and serum were spiked with different amounts of progesterone, purified as described in the Sample Purification section, serially diluted and tested in the assay. The results are shown below. The error bars represent standard deviations obtained from multiple dilutions of each sample.

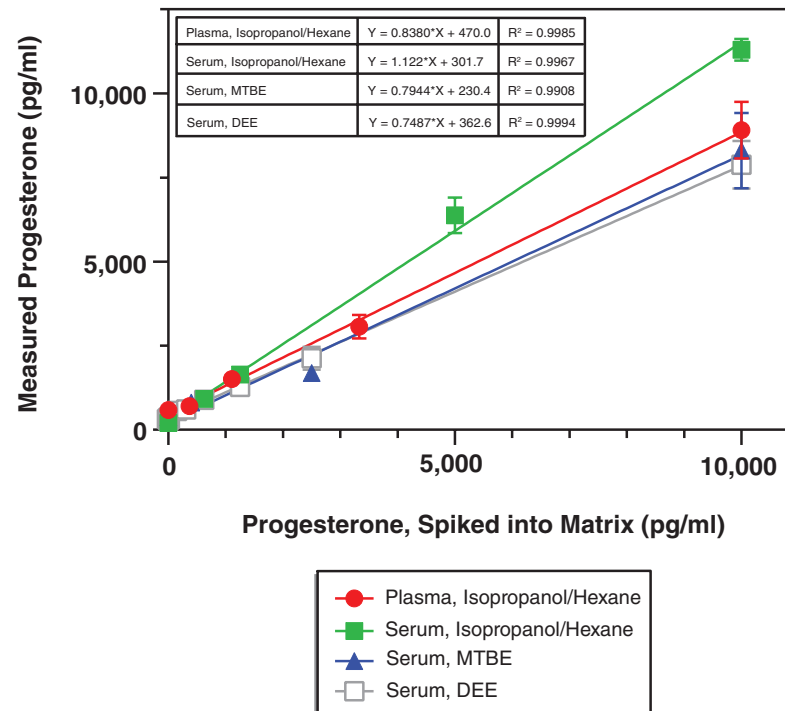


Figure 3. Spike and recovery of Progesterone in plasma and serum

Parallelism

Plasma and serum samples were purified as described in the Sample Preparation section, then serially diluted and validated in the assay.

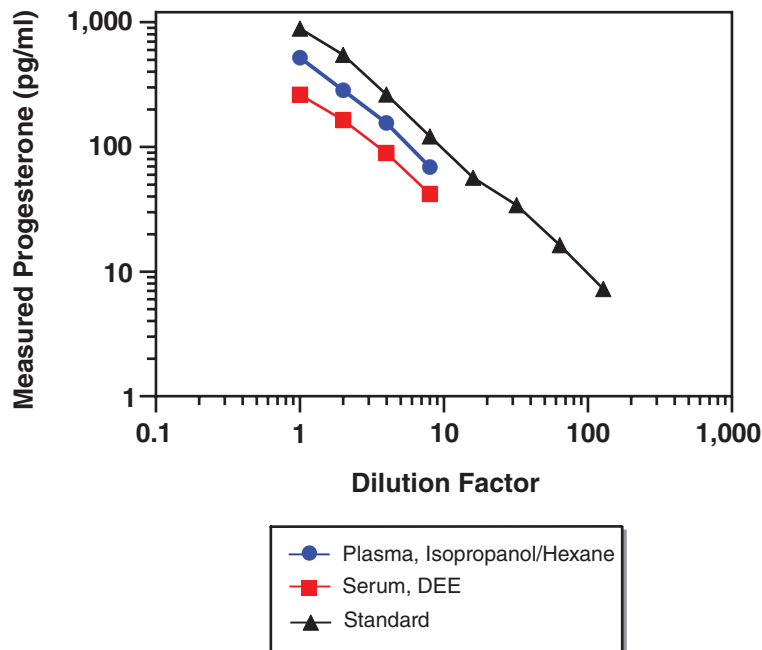


Figure 4. Parallelism in the Progesterone ELISA Kit

ASSAY PROTOCOL

Preparation of Assay-Specific Reagents

Progesterone ELISA Standard

Equilibrate a pipette tip in ethanol by repeatedly filling and expelling the tip with ethanol several times. Using the equilibrated pipette tip, transfer 100 μ l of the Progesterone ELISA Standard (Item No. 482604) into a clean test tube, then dilute with 900 μ l UltraPure water. The concentration of this solution (the bulk standard) will be 10 ng/ml.

NOTE: If assaying culture medium samples that have not been diluted with ELISA Buffer, culture medium should be used in place of ELISA Buffer for dilution of the standard curve.

To prepare the standard for use in ELISA: Obtain eight clean test tubes and number them #1 through #8. Aliquot 900 μ l ELISA Buffer to tube #1 and 500 μ l ELISA Buffer to tubes #2-8. Transfer 100 μ l of the bulk standard (10 ng/ml) to tube #1 and mix thoroughly. The concentration of this standard, the first point on the standard curve, will be 1 ng/ml (1,000 pg/ml). Serially dilute the standard by removing 500 μ l from tube #1 and placing in tube #2; mix thoroughly. Next, remove 500 μ l from tube #2 and place it into tube #3; mix thoroughly. Repeat this process for tubes #4-8. These diluted standards should not be stored for more than 24 hours.

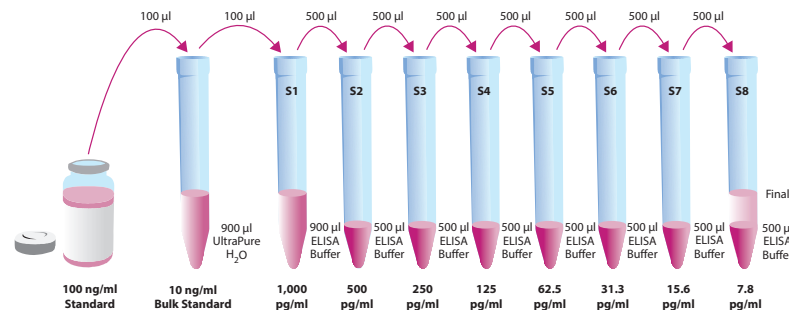


Figure 5. Preparation of the Progesterone standards

Progesterone AChE Tracer

Reconstitute the Progesterone AChE Tracer as follows:

100 dtn Progesterone AChE Tracer (96-well kit; Item No. 482600):
Reconstitute with 6 ml ELISA Buffer.

OR

500 dtn Progesterone AChE Tracer (480-well kit; Item No. 482600):
Reconstitute with 30 ml ELISA Buffer.

Store the reconstituted Progesterone AChE Tracer at 4°C (*do not freeze!*) and use within four weeks. A 20% surplus of tracer has been included to account for any incidental losses.

Progesterone ELISA Antiserum

Reconstitute the Progesterone ELISA Antiserum as follows:

100 dtn Progesterone ELISA Antiserum (96-well kit; Item No. 482602):
Reconstitute with 6 ml ELISA Buffer.

OR

500 dtn Progesterone ELISA Antiserum (480-well kit; Item No. 482602):
Reconstitute with 30 ml ELISA Buffer.

Store the reconstituted Progesterone ELISA Antiserum at 4°C. It will be stable for at least four weeks. A 20% surplus of antiserum has been included to account for any incidental losses.

Plate Set Up

The 96-well plate(s) included with this kit is supplied ready to use. It is not necessary to rinse the plate(s) prior to adding the reagents. *NOTE: If you do not need to use all the strips at once, place the unused strips back in the plate packet and store at 4°C. Be sure the packet is sealed with the desiccant inside.*

Each plate or set of strips must contain a minimum of two blanks (Blk), two non-specific binding wells (NSB), two maximum binding wells (B_0), and an eight point standard curve run in duplicate. *NOTE: Each assay must contain this minimum configuration in order to ensure accurate and reproducible results.* Each sample should be assayed at two dilutions and each dilution should be assayed in duplicate. For statistical purposes, we recommend assaying samples in triplicate.

A suggested plate format is shown in Figure 6, below. The user may vary the location and type of wells present as necessary for each particular experiment. The plate format provided below has been designed to allow for easy data analysis using a convenient spreadsheet offered by Cayman (see page 23, for more details). We suggest you record the contents of each well on the template sheet provided (see page 33).

	1	2	3	4	5	6	7	8	9	10	11	12
A	Blk	S1	S1	1	1	1	9	9	9	17	17	17
B	Blk	S2	S2	2	2	2	10	10	10	18	18	18
C	NSB	S3	S3	3	3	3	11	11	11	19	19	19
D	NSB	S4	S4	4	4	4	12	12	12	20	20	20
E	B_0	S5	S5	5	5	5	13	13	13	21	21	21
F	B_0	S6	S6	6	6	6	14	14	14	22	22	22
G	B_0	S7	S7	7	7	7	15	15	15	23	23	23
H	TA	S8	S8	8	8	8	16	16	16	24	24	24

Blk - Blank
TA - Total Activity
NSB - Non-Specific Binding
 B_0 - Maximum Binding
S1-S8 - Standards 1-8
1-24 - Samples

Figure 6. Sample plate format

Performing the Assay

Pipetting Hints
<ul style="list-style-type: none">Use different tips to pipette each reagent.Before pipetting each reagent, equilibrate the pipette tip in that reagent (i.e., slowly fill the tip and gently expel the contents, repeat several times).Do not expose the pipette tip to the reagent(s) already in the well.

Addition of the Reagents

1. ELISA Buffer

Add 100 µl ELISA Buffer to NSB wells. Add 50 µl ELISA Buffer to B₀ wells. If culture medium was used to dilute the standard curve, substitute 50 µl of culture medium for ELISA Buffer in the NSB and B₀ wells (i.e., add 50 µl culture medium to NSB and B₀ wells and 50 µl ELISA Buffer to NSB wells).

2. Progesterone ELISA Standard

Add 50 µl from tube #8 to both of the lowest standard wells (S8). Add 50 µl from tube #7 to each of the next two standard wells (S7). Continue with this procedure until all the standards are aliquoted. The same pipette tip should be used to aliquot all the standards. Before pipetting each standard, be sure to equilibrate the pipette tip in that standard.

3. Samples

Add 50 µl of sample per well. Each sample should be assayed at a minimum of two dilutions. Each dilution should be assayed in duplicate (triplicate recommended).

4. Progesterone AChE Tracer

Add 50 µl to each well except the TA and the Blk wells.

5. Progesterone ELISA Antiserum

Add 50 µl to each well except the TA, the NSB, and the Blk wells.

Well	ELISA Buffer	Standard/Sample	Tracer	Antiserum
Blk	-	-	-	-
TA	-	-	5 µl (at devl. step)	-
NSB	100 µl	-	50 µl	-
B ₀	50 µl	-	50 µl	50 µl
Std/Sample	-	50 µl	50 µl	50 µl

Table 1. Pipetting summary

Incubation of the Plate

Cover each plate with plastic film (Item No. 400012) and incubate for one hour at room temperature on an orbital shaker.

Development of the Plate

- Reconstitute Ellman’s Reagent immediately before use (20 ml of reagent is sufficient to develop 100 wells):

100 dtn vial Ellman’s Reagent (96-well kit; Item No. 400050): Reconstitute with 20 ml of UltraPure water.

OR

250 dtn vial Ellman’s Reagent (480-well kit; Item No. 400050): Reconstitute with 50 ml of UltraPure water.

NOTE: Reconstituted Ellman’s Reagent is unstable and should be used the same day it is prepared; protect the Ellman’s Reagent from light when not in use. Extra vials of the reagent have been provided should a plate need to be re-developed or multiple assays run on different days.

2. Empty the wells and rinse five times with Wash Buffer.
3. Add 200 μ l of Ellman's Reagent to each well.
4. Add 5 μ l of tracer to the TA wells.
5. Cover the plate with plastic film. Optimum development is obtained by using an orbital shaker equipped with a large, flat cover to allow the plate(s) to develop in the dark. This assay typically develops (i.e., B_0 wells ≥ 0.3 A.U. (blank subtracted)) in 60-90 minutes.

Reading the Plate

1. Wipe the bottom of the plate with a clean tissue to remove fingerprints, dirt, etc.
2. Remove the plate cover being careful to keep Ellman's Reagent from splashing on the cover. *NOTE: Any loss of Ellman's Reagent will affect the absorbance readings. If Ellman's Reagent is present on the cover, use a pipette to transfer the Ellman's Reagent into the well. If too much Ellman's Reagent has splashed on the cover to easily redistribute back into the wells, wash the plate three times with wash buffer and repeat the development with fresh Ellman's Reagent.*
3. Read the plate at a wavelength between 405 and 420 nm. The absorbance may be checked periodically until the B_0 wells have reached a minimum of 0.3 A.U. (blank subtracted). The plate should be read when the absorbance of the B_0 wells are in the range of 0.3-1.5 A.U. (blank subtracted). If the absorbance of the wells exceeds 2.0, wash the plate, add fresh Ellman's Reagent and let it develop again.

ANALYSIS

Many plate readers come with data reduction software that plot data automatically. Alternatively a spreadsheet program can be used. The data should be plotted as either $\%B/B_0$ versus log concentration using a four-parameter logistic fit or as logit B/B_0 versus log concentration using a linear fit. *NOTE: Cayman has a computer spreadsheet available for data analysis. Please contact Technical Service or visit our website (www.caymanchem.com/analysis/elisa) to obtain a free copy of this convenient data analysis tool.*

Calculations

Preparation of the Data

The following procedure is recommended for preparation of the data prior to graphical analysis.

NOTE: If the plate reader has not subtracted the absorbance readings of the blank wells from the absorbance readings of the rest of the plate, be sure to do that now.

1. Average the absorbance readings from the NSB wells.
2. Average the absorbance readings from the B_0 wells.
3. Subtract the NSB average from the B_0 average. This is the corrected B_0 or corrected maximum binding.
4. Calculate the B/B_0 (Sample or Standard Bound/Maximum Bound) for the remaining wells. To do this, subtract the average NSB absorbance from the S1 absorbance and divide by the corrected B_0 (from Step 3). Repeat for S2-S8 and all sample wells. (To obtain $\%B/B_0$ for a logistic four-parameter fit, multiply these values by 100.)

NOTE: The TA values are not used in the standard curve calculations. Rather, they are used as a diagnostic tool; the corrected B_0 divided by the actual TA (10X measured absorbance) will give the %Bound. This value should approximate the %Bound that can be calculated from the Sample Data (see page 26). Erratic absorbance values and a low (or no) %Bound could indicate the presence of organic solvents in the buffer or other technical problems (see page 30 for Troubleshooting).

Plot the Standard Curve

Plot %B/B₀ for standards S1-S8 *versus* Progesterone concentration using linear (y) and log (x) axes and perform a 4-parameter logistic fit.

Alternative Plot - The data can also be linearized using a logit transformation. The equation for this conversion is shown below. *NOTE: Do not use %B/B₀ in this calculation.*

$$\text{logit (B/B}_0\text{)} = \ln [\text{B/B}_0\text{}/(1 - \text{B/B}_0\text{])}$$

Plot the data as logit (B/B₀) *versus* log concentrations and perform a linear regression fit.

Determine the Sample Concentration

Calculate the B/B₀ (or %B/B₀) value for each sample. Determine the concentration of each sample using the equation obtained from the standard curve plot. *NOTE: Remember to account for any concentration or dilution of the sample prior to the addition to the well.* Samples with %B/B₀ values greater than 80% or less than 20% should be re-assayed as they generally fall out of the linear range of the standard curve. A 20% or greater disparity between the apparent concentration of two different dilutions of the same sample indicates interference which could be eliminated by purification.

Performance Characteristics

Sample Data

The standard curve presented here is an example of the data typically produced with this kit; however, your results will not be identical to these. You **must** run a new standard curve. Do not use the data below to determine the values of your samples. Your results could differ substantially.

	Raw Data		Average	Corrected
Total Activity	0.701	0.703	0.702	
NSB	-0.001	-0.002	-0.002	
B ₀	0.967	0.985		
	0.971	0.964	0.972	0.974

Dose (pg/ml)	Raw Data		Corrected		%B/B ₀	
1,000	0.135	0.136	0.137	0.138	14.1	14.1
500	0.196	0.202	0.198	0.204	20.3	20.9
250	0.279	0.279	0.281	0.281	28.8	28.8
125	0.390	0.389	0.392	0.391	40.2	40.1
62.5	0.497	0.510	0.499	0.512	51.2	52.5
31.3	0.609	0.606	0.611	0.608	62.8	62.4
15.6	0.723	0.717	0.725	0.719	74.5	73.9
7.8	0.783	0.796	0.785	0.798	80.6	82.0

Table 2. Typical results

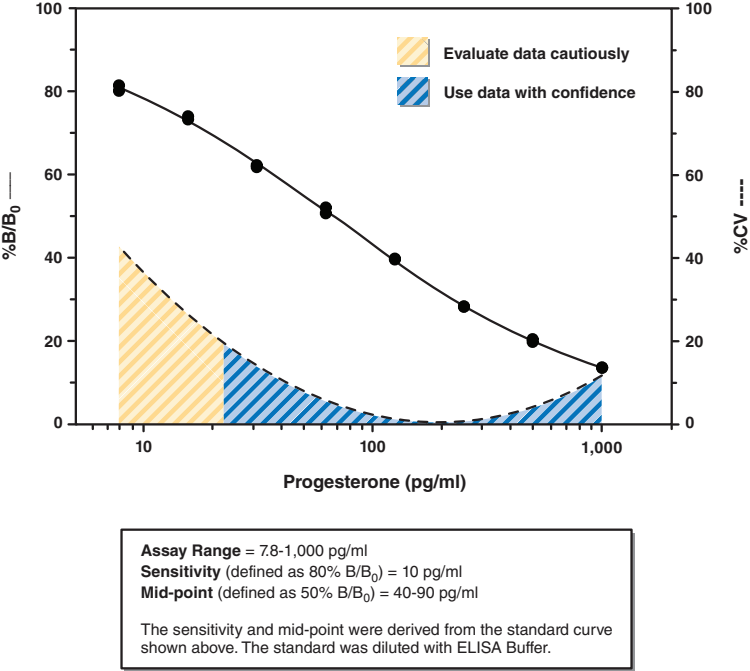


Figure 7. Typical standard curve

Precision:

The intra- and inter-assay CVs have been determined at multiple points on the standard curve. These data are summarized in the graph on page 27 and in the table below.

Dose (pg/ml)	%CV*	
	Intra-assay variation	Inter-assay variation
1,000	7.3	16.4
500	6.9	13.8
250	4.9	1.5
125	7.5	2.9
62.5	7.5	12.5
31.3	10.7	7.7
15.6	12.1	14.3
7.8	54.5	7.7

Table 3. Intra- and inter-assay variation

*%CV represents the variation in concentration (not absorbance) as determined using a reference standard curve.

Cross Reactivity:

Compound	Cross Reactivity
Progesterone	100%
Pregnenolone	14.0%
17β-Estradiol	7.2%
5β-Pregnan-3α-ol-20-one	6.7%
17α-Hydroxyprogesterone	3.6%
17α-Hydroxypregnenolone	0.44%
Corticosterone	0.26%
Testosterone	<0.05%
17α-Estradiol	<0.01%
Estriol	<0.01%
Estrone	<0.01%
5α-Pregnan-3α,20α-diol	<0.01%
5β-Pregnan-3α,20α-diol	<0.01%
5β-Pregnan-3α,20α-diol glucuronide	<0.01%
Promegestone	<0.01%
DHEA	0.015%
Cortisol	0.006%
DHEA Sulfate	0.006%

Table 4. Cross Reactivity of the Progesterone ELISA

RESOURCES

Troubleshooting

Problem	Possible Causes	Recommended Solutions
Erratic values; dispersion of duplicates	A. Trace organic contaminants in the water source B. Poor pipetting/technique	A. Replace activated carbon filter or change source of UltraPure water
High NSB (>10% of B ₀)	A. Poor washing B. Exposure of NSB wells to specific antibody	Rewash plate and redevelop
Very low B ₀	A. Trace organic contaminants in the water source B. Plate requires additional development time C. Dilution error in preparing reagents	A. Replace activated carbon filter or change source of UltraPure water B. Return plate to shaker and re-read later
Low sensitivity (shift in dose response curve)	Standard is degraded	Replace standard
Analyses of two dilutions of a biological sample do not agree (i.e., more than 20% difference)	Interfering substances are present	Purify sample prior to analysis by ELISA ⁵
Only Total Activity (TA) wells develop	Trace organic contaminants in the water source	Replace activated carbon filter or change source of UltraPure water

Additional Reading

Go to www.caymanchem.com/582601/references for a list of publications citing the use of Cayman Chemical's PGE₂ ELISA Kit.

References

1. Erickson, G.F. The ovary: Basic principles and concepts. A. Physiology, Chapter 17, in Endocrinology and metabolism. Felig, P., Baxter, J.D., and Frohman, L.A., editors, 3rd ed., McGraw-Hill, Inc. Health Professions Division, New York, 973-1015 (1995).
2. Tallon, D.F., Gosling, J.P., Buckley, P.M., *et al.* Direct solid-phase enzyme immunoassay of progesterone in saliva. *Clin. Chem.* **30**, 1507-1511 (1984).

3. Maclouf, J., Grassi, J., and Pradelles, P. Development of enzyme-immunoassay techniques for the measurement of eicosanoids, Chapter 5, *in* Prostaglandin and Lipid Metabolism in Radiation Injury. Walden, T.L., Jr. and Hughes, H.N., editors, Plenum Press, Rockville, 355-364 (1987).

4. Pradelles, P., Grassi, J., and Maclouf, J.A. Enzyme immunoassays of eicosanoids using acetylcholinesterase as label: An alternative to radioimmunoassay. *Anal. Chem.* **57**, 1170-1173 (1985).

5. Maxey, K.M., Maddipati, K.R., and Birkmeier, J. Interference in enzyme immunoassays. *J. Clin. Immunoassay* **15**, 116-120 (1992).

12								
11								
10								
9								
8								
7								
6								
5								
4								
3								
2								
1								
	A	B	C	D	E	F	G	H

Warranty and Limitation of Remedy

Buyer agrees to purchase the material subject to Cayman's Terms and Conditions. Complete Terms and Conditions including Warranty and Limitation of Liability information can be found on our website.

This document is copyrighted. All rights are reserved. This document may not, in whole or part, be copied, photocopied, reproduced, translated, or reduced to any electronic medium or machine-readable form without prior consent, in writing, from Cayman Chemical Company.

©11/05/2020, Cayman Chemical Company, Ann Arbor, MI, All rights reserved.
Printed in U.S.A.

