SZABO SCANDIC

Part of Europa Biosite

Produktinformation

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mail@szabo-scandic.com

www.szabo-scandic.com

[linkedin.com/company/szaboscandic](https://www.linkedin.com/company/szaboscandic) 

Technically
Speaking

CEDARLANE[®] 
www.cedarlanelabs.com

Conveniently Delivering You Today's Innovations
for the Science of Tomorrow™

**LOW-TOX[®] GUINEA PIG COMPLEMENT
(For Use with Mouse Lymphocytes)**

CL4051

DESCRIPTION:

Cedarlane Low-Tox[®] Guinea Pig Complement is a highly potent source of guinea pig complement, which has been absorbed to remove background toxicity to mouse lymphocytes. Each batch is tested for high complement activity and low background toxicity to mouse thymocytes. For optimal results, use Low-Tox[®] Guinea Pig Complement at a minimum dilution of 1:20.

PREPARATION: 1 ml, lyophilized

STORAGE: Store at -20°C before reconstitution. After reconstitution, store at 4°C for a maximum of **one hour**.

RECONSTITUTION:

Reconstitute with 1 ml of ice cold distilled water. Use within **one hour** of reconstitution, keeping cold until used. If this material is refrozen, there may be a reduction in potency. If refrozen, store at -70°C.

NOTE: The reconstituted material will appear cloudy. This is a result of the lyophilization process and does not affect activity.

STERILITY:

This reagent is not sold as sterile, but can be sterilized by filtration if necessary. Dilute to working dilution and then filter to avoid loss of product.

**FOR RESEARCH USE ONLY
NOT FOR HUMAN DRUG OR DIAGNOSTIC USE**

[®] Registered Trademark of Cedarlane

rev/lh/nov/98

Visit our website for your local distributor.

CEDARLANE[®] 

www.cedarlanelabs.com

An ISO 9001:2000 and ISO 13485:2003
registered company.

In CANADA: **Toll Free: 1-800-268-5058**

4410 Paletta Court, Burlington, ON L7L 5R2 ph: (289) 288-0001, fax: (289) 288-0020
e-mail: general@cedarlanelabs.com

In the USA: **Toll Free: 1-800-721-1644**

1210 Turrentine Street, Burlington, NC 27215 ph: (336) 513-5135, fax: (336) 513-5138
e-mail: service@cedarlanelabs.com