SZABO SCANDIC

Part of Europa Biosite

Produktinformation

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mail@szabo-scandic.com

www.szabo-scandic.com

[linkedin.com/company/szaboscandic](https://www.linkedin.com/company/szaboscandic) 

TECHNICALLY Speaking

Place your order with CEDARLANE® or your local distributor.

Please contact CEDARLANE® for lot specific information.

LOW-TOX® GUINEA PIG COMPLEMENT (For Use with Mouse Lymphocytes)

CL4055

DESCRIPTION:

Cedarlane Low-Tox® Guinea Pig Complement is a highly potent source of guinea pig complement that has been absorbed with agarose to remove background toxicity to mouse lymphocytes. Each batch is tested for high complement activity and low background toxicity to mouse thymocytes. *For optimal results, use Low-Tox® Guinea Pig Complement at a minimum dilution of 1:20.*

PREPARATION: 5 ml, lyophilized

STORAGE: Store at -20°C or below before reconstitution. After reconstitution, store at 4°C for a maximum of one hour.

RECONSTITUTION:

Reconstitute with 5 ml of ice cold distilled water. Use within **one hour** of reconstitution, keeping cold until used. If this material is refrozen, there may be a reduction in potency. If refrozen, store at -70°C.

NOTE: The reconstituted material will appear cloudy. This is a result of the lyophilization process and does not affect activity.

STERILITY:

This reagent is not sold as sterile, but can be sterilized by filtration if necessary. Dilute to working dilution and then filter to avoid loss of product.

FOR LABORATORY RESEARCH ONLY

NOT FOR HUMAN DRUG OR DIAGNOSTIC USE

® Registered Trademark of Cedarlane Laboratories Ltd.

Rev/da/jun/00

For more information or to place an order please contact...

CEDARLANE®
LABORATORIES LIMITED

toll free: 1-800-268-5058
in North America

phone: (905) 878-8891 • fax: (905) 878-7800

5516 - 8th Line, R.R.#2, Hornby, Ontario, CANADA L0P 1E0

or visit our website for a list of our international distributors including contact information

website: www.cedarlanelabs.com • e-mail: info@cedarlanelabs.com