SZABO SCANDIC

Part of Europa Biosite

Produktinformation

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mail@szabo-scandic.com

www.szabo-scandic.com

[linkedin.com/company/szaboscandic](https://www.linkedin.com/company/szaboscandic) 

Technically
Speaking

CEDARLANE[®] 
www.cedarlanelabs.com

Conveniently Delivering You Today's Innovations
for the Science of Tomorrow™

**Human COMPLEMENT
(Pooled Normal Human Serum)**

CL6600-1
CL6600-10
CL6600-50
CL6600-100

DESCRIPTION:

Cedarlane's Human Complement is aseptically collected from healthy individuals.

PRESENTATION:

CL6600-1:	1 ml, frozen
CL6600-10:	10 ml, frozen
CL6600-50:	50 ml, frozen
CL6600-100:	100 ml, frozen

STORAGE AND USE:

Store at -70°C. Thaw the complement by incubating in an **ice bath**. Do not subject complement to temperatures above 4°C.. Use within one hour of thawing. If complement is to be used and refrozen, aliquot immediately into single use quantities and freeze quickly at -70°C. Store at -70°C until use. Do not refreeze. Aliquots should be dated and the activity checked before use.

COMMENTS:

All materials tested and found to be non-reactive for HBsAG, HIV, HIV-1 p24 antigen and HCV at the time of donation or thereafter by currently approved U.S.A. FDA methods. All blood products should be treated as potentially infectious source material from which this product was derived. No known test method can offer assurance that products derived from human blood will not transmit infectious agents.

Product is to be used by laboratory-trained staff. Disposal is in accordance with local regulations.

FOR RESEARCH USE ONLY

Visit our website for your local distributor.

CEDARLANE[®] 

www.cedarlanelabs.com

An ISO 9001:2000 and ISO 13485:2003
registered company.

In CANADA: **Toll Free: 1-800-268-5058**

4410 Paletta Court, Burlington, ON L7L 5R2 ph: (289) 288-0001, fax: (289) 288-0020
e-mail: general@cedarlanelabs.com

In the USA: **Toll Free: 1-800-721-1644**

1210 Turrentine Street, Burlington, NC 27215 ph: (336) 513-5135, fax: (336) 513-5138
e-mail: service@cedarlanelabs.com