# SZABO SCANDIC

Part of Europa Biosite

## Produktinformation

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[mail@szabo-scandic.com](mailto:mail@szabo-scandic.com)

[www.szabo-scandic.com](http://www.szabo-scandic.com)

[linkedin.com/company/szaboscandic](https://www.linkedin.com/company/szaboscandic) 

Technically  
Speaking

**CEDARLANE**<sup>®</sup>   
[www.cedarlanelabs.com](http://www.cedarlanelabs.com)

Conveniently Delivering You Today's Innovations  
for the Science of Tomorrow™

## Solution Pepsin

**CLSG80108**  
**Lot: W412**

**Intended Use:** This product is designed to unmask immunoreactive sites altered by either fixation or embedding media.

**Introduction:** Fixation is one of the most critical aspects of immunostaining. If the antigen is not properly fixed it will be washed out of the specimen. On the other hand, over fixation can cause severe problems such as masking or denaturation of the antigen. Formaldehyde is probably the best all-around fixative, due to its cross-linking characteristics. Length of fixation is very critical to prevent antigen masking. Immunoglobulins are especially susceptible to over fixation with formalin. In such cases, treatment with proteolytic enzyme is required to digest excess aldehyde linkages and to expose the antigen.

**Format:** 50 mL clear ready-to-use solution

**Storage:** Store at 2-8°C. Do not use beyond the expiration date stated on the label.

**Digestion Procedure:**

1. Deparaffinize tissue sections and bring them to wash buffer.
2. Remove excess buffer from slide(s) without letting the tissue dry.
3. Depending on the size of the tissue section add 1-2 drops of pepsin solution.
4. Incubate slide(s) for 5-10 minutes at 37°C.
5. Wash slide(s) with wash buffer and proceed with immunostaining.

BA 02/19/13

**Laboratory Reagent For Research Use Only**

Visit our website for your local distributor.

**CEDARLANE**<sup>®</sup> 

[www.cedarlanelabs.com](http://www.cedarlanelabs.com)

An ISO 9001:2000 and ISO 13485:2003  
registered company.

**In CANADA: Toll Free: 1-800-268-5058**

4410 Paletta Court, Burlington, ON L7L 5R2 ph: (289) 288-0001, fax: (289) 288-0020  
e-mail: [general@cedarlanelabs.com](mailto:general@cedarlanelabs.com)

**In the USA: Toll Free: 1-800-721-1644**

1210 Turrentine Street, Burlington, NC 27215 ph: (336) 513-5135, fax: (336) 513-5138  
e-mail: [service@cedarlanelabs.com](mailto:service@cedarlanelabs.com)