# SZABO SCANDIC

Part of Europa Biosite

## Produktinformation

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[mail@szabo-scandic.com](mailto:mail@szabo-scandic.com)

[www.szabo-scandic.com](http://www.szabo-scandic.com)

[linkedin.com/company/szaboscandic](https://www.linkedin.com/company/szaboscandic) 

# TECHNICALLY *Speaking*

Place your order with CEDARLANE® or your local distributor.  
Please contact CEDARLANE® for lot specific information.

## Peroxidase Block

**CLSG80120**

**Lot:**

**Size:** 50 ml

**Intended Use:** This product is designed for use in peroxidase based immunohistochemical staining procedures done on paraffin tissue sections to suppress non-specific staining due to endogenous peroxidase activity.

**Introduction:** Endogenous peroxidase activity has been observed in tissue sections during immunoperoxidase staining. To carry out immunoperoxidase staining, endogenous peroxidase must be suppressed from the tissue sections. If not properly suppressed, endogenous peroxidase can produce false immunostaining signals as well as can create high background during the immunostaining. Treatment of paraffin tissue sections with peroxidase block prior to the application of primary antibody will suppress endogenous peroxidase activity.

**Storage:** Store refrigerated. Do not freeze.

**Procedure:** Incubate paraffin embedded tissue sections for 5-10 minutes at room temperature prior to applying primary antibody. Rinse tissue with wash buffer after the incubation with peroxidase block and proceed to the next step.

**Laboratory Reagent For Research Use Only**

JV 11/01/07

For more information or to place an order please contact...

**CEDARLANE®**  
**LABORATORIES LIMITED**

*toll free: 1-800-268-5058*  
*in North America*

phone: (905) 878-8891 • fax: (905) 878-7800

5516 - 8th Line, R.R.#2, Hornby, Ontario, CANADA L0P 1E0

or visit our website for a list of our international distributors including contact information  
website: [www.cedarlanelabs.com](http://www.cedarlanelabs.com) • e-mail: [info@cedarlanelabs.com](mailto:info@cedarlanelabs.com)