



SZABO SCANDIC

Part of Europa Biosite

Produktinformation



Forschungsprodukte & Biochemikalien



Zellkultur & Verbrauchsmaterial



Diagnostik & molekulare Diagnostik



Laborgeräte & Service

Weitere Information auf den folgenden Seiten!
See the following pages for more information!



Lieferung & Zahlungsart

siehe unsere [Liefer- und Versandbedingungen](#)

Zuschläge

- Mindermengenzuschlag
- Trockeneiszuschlag
- Gefahrgutzuschlag
- Expressversand

SZABO-SCANDIC HandelsgmbH

Quellenstraße 110, A-1100 Wien

T. +43(0)1 489 3961-0

F. +43(0)1 489 3961-7

mail@szabo-scandic.com

www.szabo-scandic.com

[linkedin.com/company/szaboscandic](https://www.linkedin.com/company/szaboscandic) 

Technically
Speaking

CEDARLANE[®]
www.cedarlanelabs.com



Conveniently Delivering You Today's Innovations
for the Science of Tomorrow™

Fontana-Masson Stain Kit

CLSG80154

Lot:

Kit Components Included:

125 ml	0.2% Gold Chloride Sol'n	2-8°C
5 x 9 ml	10% Silver Nitrate Sol'n	2-8°C
125 ml	5% Sodium Thiosulfate Sol'n	Room Temp.
125 ml	Nuclear Fast Red Sol'n	Room Temp.

Description: The Fontana-Masson Stain Kit is intended for use in the histological visualization of Argentaffin cells and Melanin in paraffin or frozen sections.

Control Tissue: Any paraffin embedded tissue that contains Hair Follicles or Skin for Melanin. Small Intestine for Argentaffin.

Precautions: Avoid contact with skin and eyes.
Harmful if swallowed.
Follow all Federal, State, and local regulations regarding disposal.
Use in chemical fume hood whenever possible.

Preparation of Reagent Prior to Beginning:

Prepare Ammoniacal Silver Solution immediately prior to use.

In new or chemically cleaned glassware, mix 27ml Distilled/Deionized water with one vial of Silver Nitrate Solution (10%) and blend completely. Carefully add Concentrated Ammonium Hydroxide (Not included) one drop at a time, swirling gently after each drop. Initially the mixture will turn dark brown and then gradually become transparent with a fine layer of sediment. The solution is ready for immediate use when all sediment dissolves.

Continued...

Visit our website for your local distributor.

CEDARLANE[®]



www.cedarlanelabs.com

An ISO 9001:2000 and ISO 13485:2003
registered company.

In CANADA: Toll Free: 1-800-268-5058

4410 Paletta Court, Burlington, ON L7L 5R2 ph: (289) 288-0001, fax: (289) 288-0020
e-mail: general@cedarlanelabs.com

In the USA: Toll Free: 1-800-721-1644

1210 Turrentine Street, Burlington, NC 27215 ph: (336) 513-5135, fax: (336) 513-5138
e-mail: service@cedarlanelabs.com

Procedure:

1. Deparaffinize sections if necessary and hydrate to distilled water.
2. Place freshly mixed Ammoniacal Silver Solution in a 58-60°C waterbath and allow adequate time for temperature to equilibrate.
3. Incubate slide in warmed Ammoniacal Silver Solution for 30-60 minutes or until tissue section becomes yellowish/brown in color. (NOTE: Melanin typically stains in 30 minutes while Argentaffin stains in 50-60 minutes)
4. Rinse in several changes of distilled water.
5. Incubate slide in Gold Chloride Solution (0.2%) for 30 seconds.
6. Rinse in several changes of distilled water.
7. Incubate slide in Sodium Thiosulfate Solution (5%) for 1-2 minutes.
8. Rinse for 2 minutes in running tap water followed by 2 changes of distilled water.
9. Incubate slide in Nuclear Fast Red Solution for 5 minutes.
10. Rinse for 2 minutes in running tap water followed by 2 changes of distilled water.
11. Dehydrate very quickly in 3 changes of fresh Absolute Alcohol.
12. Clear, and mount in synthetic resin.

Reference: Sheenan, D.C., Hrapchak, B.B. Theory and Practice of Histotechnology, 2nd Edition. Battelle Press, Columbus, OH.

For Research Use Only. Not for Diagnostic or Therapeutic Work.