



SZABO SCANDIC

Part of Europa Biosite

Produktinformation



Forschungsprodukte & Biochemikalien



Zellkultur & Verbrauchsmaterial



Diagnostik & molekulare Diagnostik



Laborgeräte & Service

Weitere Information auf den folgenden Seiten!
See the following pages for more information!



Lieferung & Zahlungsart

siehe unsere [Liefer- und Versandbedingungen](#)

Zuschläge

- Mindermengenzuschlag
- Trockeneiszuschlag
- Gefahrgutzuschlag
- Expressversand

SZABO-SCANDIC HandelsgmbH

Quellenstraße 110, A-1100 Wien

T. +43(0)1 489 3961-0

F. +43(0)1 489 3961-7

mail@szabo-scandic.com

www.szabo-scandic.com

[linkedin.com/company/szaboscandic](https://www.linkedin.com/company/szaboscandic) 

SARS-CoV-2 Pseudoviral Particles, Omicron-BJ.1 Sub-variant

CATALOG NUMBER: SCV2-PsV-BJ1, 5 mL

Description

It has been known that SARS coronavirus 2 (SARS-CoV-2) uses human ACE2 as entry receptor and human proteases as entry activators. The virus surface spike protein (S) mediates SARS-CoV-2 entry into cells. To fulfill its function, SARS-CoV-2 spike binds to its human ACE2 (hACE2) receptor through its receptor-binding domain (RBD) and is proteolytically activated by human proteases.

Our **SARS-CoV-2 Omicron BJ.1 Pseudoviral Particles** are replication-deficient MLV pseudotyped with the SARS-CoV-2 spike protein of the Omicron BJ.1 sub-variant (Genbank sequence ID: UXU98303.1). They also contain the ORF for firefly luciferase as a reporter. They establish a pseudovirus cell entry assay mediated by the SARS-CoV-2 spike protein that can be conveniently measured via luciferase reporter activity. This pseudovirus assay isolates the SARS-CoV-2 viral entry from other steps of the viral infection cycle.

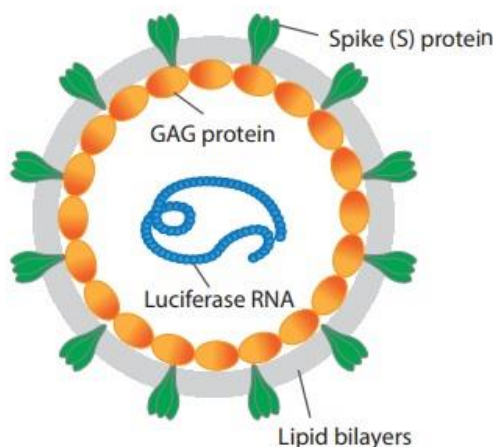


Figure 1. Illustration of the replication-deficient MLV particle pseudotyped with SARS-CoV-2 Spike protein

Applications

Our Pseudovirus Particles (PP) generate robust chemiluminescent signals in cell assays when coupled with our firefly luciferase assay kit (Catalog # [CA-L165](#)), useful for: 1) screening potential inhibitor to block SARS-CoV-2 entry and viral protein translation; 2) measuring the activity of and screening for neutralizing antibody against SARS-CoV-2-Omicron BFJ.1 (refer to [the Neutralization Assay Application Note](#)).

Features

- **Robust:** Excellent signal to noise (basal) ratio
- **Easy to use:** Amenable to HTS format (96-well, 384-well and 1536-well format)

Contents

5 ml, for one multi-well plate; PP per mL > 1.0E+07

Storage

Upon receiving it, store at -70 °C right away. Thaw* before immediate use.

***Note: read the instruction for thawing in the following protocol carefully. Do not aliquot and refreeze.**

Shelf Life:

Six months from the date of shipping when stored at -70 °C

ASSAY PROTOCOL

Note: requires a luciferase assay reagent (Catalog # [CA-L165](#)).

Cell Infection:

1. Count HEK293-ACE2 cells (Catalog # [CL-hACE2-001](#)) to be infected and seed ~20K cells per well into 96-well plates (50 μ l per well) DMEM with 10% HyClone[™] FetalClone[™] II Serum (no antibiotics) or 5K cells per well into 384-well plates (15 μ l per well).
2. Culture cells overnight to make sure the cells stably adhere to the plates.
3. On the 2nd day, remove media, add 50 μ l SARS-CoV-2 Omicron BJ.1 Pseudoviral Particles* into each well (12.5 μ l for 384-well plate). Spin at 700 rpm for 15 min at 4°C.

****Note: thaw the pseudoviral particles quickly in the room temperature water (< 30 minutes, do not shake) and use right away. Discard the unused portion (do not re-freeze or leave it on ice for later use).***

4. Incubate for 2 hrs at 37 °C.
5. Add 50 μ l DMEM with 10% FC into each well (12.5 μ l for 384-well plates).
6. Incubate for 42 hrs at 37 °C.

Measurement of Luciferase Activity in Infected cells

1. Do not remove supernatant. Add 100 μ l eEnzyme's luciferase assay reagent (25 μ l for 384-well plates).
2. Read in a luminescence plate reader and record the data.

