



# SZABO SCANDIC

Part of Europa Biosite

## Produktinformation



Forschungsprodukte & Biochemikalien



Zellkultur & Verbrauchsmaterial



Diagnostik & molekulare Diagnostik



Laborgeräte & Service

Weitere Information auf den folgenden Seiten!  
See the following pages for more information!



### Lieferung & Zahlungsart

siehe unsere [Liefer- und Versandbedingungen](#)

### Zuschläge

- Mindermengenzuschlag
- Trockeneiszuschlag
- Gefahrgutzuschlag
- Expressversand

### SZABO-SCANDIC HandelsgmbH

Quellenstraße 110, A-1100 Wien

T. +43(0)1 489 3961-0

F. +43(0)1 489 3961-7

[mail@szabo-scandic.com](mailto:mail@szabo-scandic.com)

[www.szabo-scandic.com](http://www.szabo-scandic.com)

[linkedin.com/company/szaboscandic](https://www.linkedin.com/company/szaboscandic) 

# EconoTek HRP Anti-Polyvalent (DAB) Ready-To-Use (70 slide)

**Description:** Species of Origin: Goat  
Antigen Specificity: Anti-Mouse, Rat, Rabbit, Guinea Pig  
Preadsorbed Against: Human  
Enzyme Conjugate: Peroxidase  
Chromogen Substrate: Diaminobenzidine (DAB)

**Contents:** Super Block 8 ml  
Peroxide Block 8 ml  
EconoTek Anti-Polyvalent 8 ml  
EconoTek HRP 8 ml  
DAB Chromogen 3 ml  
DAB Substrate 5 ml x 7 vials

**Uses/Limitations:** Not to be taken internally.  
For In-Vitro Diagnostic use only.  
Do not use if reagent becomes cloudy.  
Do not use past expiration date.  
Use caution when handling reagents.  
Non-Sterile.


**Storage:** Store at 2-8°C.

**Precautions:** Avoid contact with skin and eyes.  
Harmful if swallowed.  
Follow all Federal, State, and local regulations regarding disposal.

## Procedure:


1. Deparaffinize and rehydrate tissue section.
2. To reduce nonspecific background staining due to endogenous peroxidase, incubate slide in hydrogen peroxide for 5 minutes.
3. Wash 2 times in buffer.
4. If required, incubate tissue in digestive enzyme.
5. Wash 4 times in buffer.
6. Apply Super Block (blue cap), and incubate for 5-10 minutes at room temperature to block nonspecific background staining.  
*Note: Do not exceed 10 minutes or there may be a reduction in desired stain.*
7. Wash 1 time in buffer.
8. Apply primary antibody and incubate according to manufacturer's protocol.

Storage: 2° C  8° C



ScyTek Laboratories, Inc.  
205 South 600 West  
Logan, UT 84321  
U.S.A.

CE 

  
Emergo Europe  
Prinsessegracht 20  
2514 AP The Hague, The Netherlands

9. Wash 4 times in buffer.
10. Apply EconoTek Biotinylated Anti-polyvalent (yellow cap), and incubate for 30 minutes at room temperature.
11. Wash 4 times in buffer.
12. Apply EconoTek HRP (red cap), and incubate for 30 minutes at room temperature.
13. Rinse 4 times in buffer.
14. Add 8 drops of DAB chromogen to one 5ml vial of DAB Substrate. Mix well and apply to tissue for 5 minutes.
15. Rinse 1 time in Deionized water.
16. Apply chromogen/substrate mixture to tissue for another 5 minutes.
17. Rinse in Deionized water.
18. Counterstain as desired.  
**WARNING:**DAB is a suspected carcinogen. Handle with care and dispose of according to all regulations.
19. Counterstain and coverslip

### Troubleshooting Guide

#### Overstaining:

1. Concentration of the primary antibody was too high or the incubation time was too long.
2. Temperature during incubation was too high.
3. Incubation time with link antibody or streptavidin/enzyme label was too long.

#### Nonspecific Background Staining:

1. Inadequate rinsing between steps.
2. Tissue was allowed to dry with reagents on.
3. Folds in tissue trapped reagents.
4. Tissue contains endogenous peroxidase.
5. Tissue contains endogenous biotin.
6. Antigen migrated in tissue.
7. Excessive tissue adhesive on slides.
8. Inadequate blocking with protein block.


#### Weak Staining:

1. Primary antibody concentration was too low or incubation time was too short.
2. Reagents are past their expiration date.
3. Inadequate removal of wash water between steps, resulting in dilution of reagents.
4. Room temperature was excessively cool.
5. The primary antibody does not recognize an antigen that survives fixation and embedding in high enough amounts.
6. Excessive incubation with protein block (Super Block).

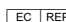
#### No Staining:

1. Steps were inadvertently left out.

Storage: 2° C  8° C


 ScyTek Laboratories, Inc.  
205 South 600 West  
Logan, UT 84321  
U.S.A.


 

  
Emergo Europe  
Prinsessegracht 20  
2514 AP The Hague, The Netherlands

P.O. Box 3286 - Logan, Utah 84323, U.S.A. - Tel. (800) 729-8350 – Tel. (435) 755-9848 - Fax (435) 755-0015 - [www.scytek.com](http://www.scytek.com)

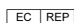
2. There is no antigen in the tissue.
3. The primary antibody is not of mouse, rat, rabbit or guinea pig origin.
4. Chromogenic substrate has been replaced with another that is not intended for use with peroxidase.

Storage: 2° C  8° C



ScyTek Laboratories, Inc.  
205 South 600 West  
Logan, UT 84321  
U.S.A.

CE 

  
Emergo Europe  
Prinsessegracht 20  
2514 AP The Hague, The Netherlands