



# SZABO SCANDIC

Part of Europa Biosite

## Produktinformation



Forschungsprodukte & Biochemikalien



Zellkultur & Verbrauchsmaterial



Diagnostik & molekulare Diagnostik



Laborgeräte & Service

Weitere Information auf den folgenden Seiten!  
See the following pages for more information!



### Lieferung & Zahlungsart

siehe unsere [Liefer- und Versandbedingungen](#)

### Zuschläge

- Mindermengenzuschlag
- Trockeneiszuschlag
- Gefahrgutzuschlag
- Expressversand

### SZABO-SCANDIC HandelsgmbH

Quellenstraße 110, A-1100 Wien

T. +43(0)1 489 3961-0

F. +43(0)1 489 3961-7

[mail@szabo-scandic.com](mailto:mail@szabo-scandic.com)

[www.szabo-scandic.com](http://www.szabo-scandic.com)

[linkedin.com/company/szaboscandic](https://www.linkedin.com/company/szaboscandic) 

# ImmPRESS® HRP Universal PLUS Polymer Kit

Peroxidase

(Horse Anti-Mouse/Rabbit IgG)



**Cat. No.** MP-7800

**Storage** Store reagents in original bottles at 2-8 °C. Do not freeze.

**Description** Instructions for Immunohistochemical staining using mouse or primary antibodies.

The ImmPRESS HRP Universal PLUS Polymer Kit is an enzymatic, non-biotin amplification system that produces crisp, highly sensitive, specific staining with low background.

The reagents in the ImmPRESS HRP Universal PLUS Polymer Kit are ready-to-use—no mixing or titering is necessary to obtain optimal staining. Dilution of the reagents or changes in suggested incubation time may affect performance. The reagents are supplied in convenient dropper bottles. (To remove the drop dispenser tip, press laterally with thumb until tip snaps off).

A number of different wash buffers can be used with the ImmPRESS reagents. One of the most common is 10 mM sodium phosphate, pH 7.5, 0.9% saline (PBS). 0.1% Tween® 20 detergent may be added to the wash buffer and is especially recommended for use with automated stainers.

**Kit Components**

Product Name	Volume
BLOXALL® Endogenous Blocking Solution	15 ml
Normal Horse Serum, 2.5%	15 ml
ImmPRESS HRP Universal Polymer Reagent (Horse Anti-Mouse/Rabbit IgG)	15 ml
ImmPACT® DAB EqV Reagent 1 (Chromogen)	7.5 ml
ImmPACT DAB EqV Reagent 2 (Diluent)	7.5 ml

The ImmPRESS Universal PLUS Polymer Kit (15 ml) will stain approximately 75-150 sections.

**Staining Procedure**

1. For paraffin sections, deparaffinize and hydrate tissue sections through xylenes or other clearing agents and graded alcohol series.

For frozen sections or cell preparations fix with acetone or an appropriate fixative for the antigen under study, if necessary.

Wash for 5 minutes in tap water.

2. If antigen unmasking is required, perform this procedure using a Vector® Antigen Unmasking Solution, Citrate-based, pH 6.0 (H-3300) or Tris-based, pH 9.0 (H-3301).



Human colon (FFPE): Stained for cytokeratin using a rabbit primary antibody and ImmPRESS HRP Universal PLUS Polymer Kit (MP-7800). No counterstain.

- If quenching of endogenous peroxidase activity is required, incubate the sections in BLOXALL Endogenous Blocking Solution for 10 minutes.
- Wash in buffer for 5 minutes.
- Incubate sections for 20 minutes with Normal Horse Serum, 2.5%.
- Tip off excess serum from sections.
- Incubate with mouse or rabbit primary antibody diluted in appropriate antibody diluent solution, such as diluted normal horse serum or BSA.
- Wash in buffer for 5 minutes.
- Incubate for 30 minutes with ImmPRESS HRP Universal Polymer Reagent.
- Wash for 2 x 5 minutes in buffer.
- Combine equal volumes of ImmPACT DAB EqV Reagent 1 with Reagent 2. Mix well.
- Incubate in ImmPACT DAB EqV working solution until desired stain intensity develops, approximately 2-10 minutes.
- Wash for 2 x 5 minutes in buffer.
- Rinse sections in tap water.
- Counterstain (optional), clear and mount.

Detailed product listings, specifications, protocols and additional information is available on our website: [vectorlabs.com](http://vectorlabs.com)

For research use only. Not intended for animal or human therapeutic or diagnostic use.

©2022 Vector Laboratories. All rights reserved.

BLOXALL, ImmPACT, ImmPRESS, and Vector are trademarks of Vector Laboratories, Inc.

Tween is a trademark of Atlas Chemical Industries.

MP-7800.LBL-00285.Rev01