



SZABO SCANDIC

Part of Europa Biosite

Produktinformation



Forschungsprodukte & Biochemikalien



Zellkultur & Verbrauchsmaterial



Diagnostik & molekulare Diagnostik



Laborgeräte & Service

Weitere Information auf den folgenden Seiten!
See the following pages for more information!



Lieferung & Zahlungsart

siehe unsere [Liefer- und Versandbedingungen](#)

Zuschläge

- Mindermengenzuschlag
- Trockeneiszuschlag
- Gefahrgutzuschlag
- Expressversand

SZABO-SCANDIC HandelsgmbH

Quellenstraße 110, A-1100 Wien

T. +43(0)1 489 3961-0

F. +43(0)1 489 3961-7

mail@szabo-scandic.com

www.szabo-scandic.com

[linkedin.com/company/szaboscandic](https://www.linkedin.com/company/szaboscandic) 

M.O.M.[®] (Mouse on Mouse) Elite[®] Immunodetection Kit

Peroxidase

Cat. No.	PK-2200
Storage	Store reagents in original bottles at 2–8 °C. We recommend that the reagents be kept in the box in which they were supplied. Do not freeze.
Description	<p>The Vector[®] M.O.M. (Mouse on Mouse) Immunodetection Kits are designed specifically to localize mouse primary antibodies on mouse tissues.</p> <p>The M.O.M. Elite Immunodetection Kit can be used with normal and genetically engineered mouse models, including transgenic, xenograft, knock out and other mutant strains.</p>

Kit Components

Product Name	Volume
M.O.M. Blocking Reagent	1 ml
M.O.M. Protein Concentrate	6 ml
M.O.M. Biotinylated Anti-Mouse IgG	0.1 ml
Reagent A (Avidin, ABC Elite)	1 ml
Reagent B (Biotinylated HRP, ABC Elite)	1 ml

The M.O.M. Elite Immunodetection Kit contains enough stock reagents to produce about 25 ml of working solution which is generally sufficient to stain approximately 250 tissue sections.

Preparation of M.O.M. Working Solutions


- M.O.M. Mouse IgG Blocking Reagent: add 2 drops (90 µl) of M.O.M. Blocking Reagent stock solution to 2.5 ml of PBS or TBS. †
- M.O.M. Diluent: add 600 µl of M.O.M. Protein Concentrate stock solution to 7.5 ml of PBS or TBS. ††
- M.O.M. Biotinylated Anti-Mouse IgG Reagent: add 10 µl of M.O.M. Biotinylated Anti-Mouse IgG stock solution to 2.5 ml of M.O.M. Diluent prepared above.
- VECTASTAIN[®] Elite[®] ABC Reagent: add 2 drops of Reagent A to 2.5 ml of PBS or TBS. Mix. Add 2 drops of Reagent B and mix. Allow VECTASTAIN Elite ABC Reagent to stand for 15–30 minutes prior to use.

† PBS: 10 mM sodium phosphate, 0.15 M NaCl, pH 7.4–7.8
TBS: 50 mM TRIS, 0.15 M NaCl, pH 7.5–7.8

†† Note: 7.5 ml of M.O.M. Diluent provides sufficient reagent for use in steps 8, 9, 11.

Staining Procedure

1. For paraffin sections, deparaffinize and hydrate tissue sections through xylenes or other clearing agents and graded alcohol series. For frozen sections or cell preparations, fix with acetone or an appropriate fixative for the antigen under study. Air dry. Rinse for 5 minutes in tap water.

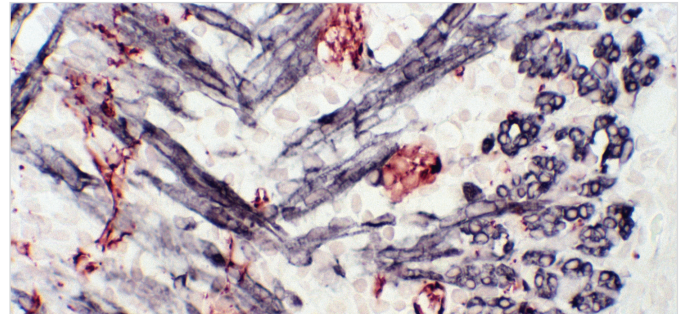

Newborn Mouse Tongue - Double Label: Synapsin (m), M.O.M. Elite[®] Immunodetection Kit, Peroxidase (PK-2200), Vector NovaRED[®] (red), Desmin (m), M.O.M. Elite[®] Immunodetection Kit, Peroxidase (PK-2200), Vector DAB-Ni (black)

2. If antigen unmasking is required, perform this procedure using an Antigen Unmasking Solution, Citrate-based, pH 6.0 (H-3300) or Tris-based, pH 9.0 (H-3301).
3. Block endogenous enzyme activity, if necessary, by incubating sections with BLOXALL[®] Blocking Solution (SP-6000) for 10 minutes.
4. Wash 2 x 2 minutes in PBS or TBS.
5. Perform Avidin/Biotin blocking if required*, using Avidin/Biotin Blocking Kit (SP-2001).
6. Incubate for 1 hour in working solution of prepared M.O.M. Mouse IgG Blocking Reagent.
7. Wash 2 x 2 minutes in PBS or TBS**.
8. Incubate for 5 minutes in working solution of prepared M.O.M. Diluent**.
9. Dilute primary antibody in M.O.M. Diluent to the appropriate concentration. Tip off excess M.O.M. Diluent from sections and apply diluted primary antibody. Incubate for 30 minutes**.
10. Wash for 2 x 2 minutes in PBS or TBS**.
11. Apply working solution of prepared M.O.M. Biotinylated Anti-Mouse IgG Reagent. Incubate sections for 10 minutes**.
12. Wash for 2 x 2 minutes in PBS or TBS.
13. Apply prepared VECTASTAIN Elite ABC Reagent. Incubate sections for 5 minutes.
14. Wash for 2 x 5 minutes in PBS or TBS.
15. Prepare and apply peroxidase substrate solution (not included in kit) according to substrate kit instructions.

* When appropriate control sections have shown that endogenous enzyme or endogenous avidin/biotin activity is not present, step 3 and/or step 5 may be omitted.

** It is recommended that the exact times described in steps 7–11 be used in the staining protocol. Longer incubation may result in an increase in background staining.

Detailed product listings, specifications, protocols and additional support information such as our Troubleshooting Guide: Mouse Antibodies on Mouse Tissue are available on our website: vectorlabs.com